# Ultrospec<sup>™</sup> 7000/8000/9000

# UV-Visible Spectrophotometers User Manual





Biochrom US 84 October Hill Rd Holliston, MA 01746-1388 USA

Telephone: 1-508-893-8999 Toll Free: 1-800-272-2775 Fax: 1-508-429-5732 support@hbiosci.com www.biochromspectros.com

# **CONTENTS**:

1.	HEALTH & SAFETY	4
		4
	1.2. Unpacking & Installation	5
	1.2.1. Instrument Connections	0
	1.3.1 Controls and Indicators	7
	1.3.2 Intended Use	7
	133 Instrument Preparation	7
	1.3.4. Post Run Procedures	8
	1.3.5. Performance Validation	8
	1.4. User Maintenance	8
	1.4.1. Troubleshooting	9
	1.5. Customer Support Contacts	10
	1.6. Service, Repair or Return	10
2		10
۷.	9000 SPECTROPHOTOMETER	11
3.	USE WITH Datrys PC SOFTWARE	12
4.	FREQUENTLY USED ICONS	13
5.	PERFORMING A MEASUREMENT	14
	5.1. Ultrospec 7000/8000/9000	14
6.	LAMP MODE	15
7.	TYPES OF BOXES	16
8	SETTINGS	17
0.	8.1. Date and time	17
	8.2. Regional	17
	8.3. Data Output	18
	8.4. User Interface	18
	8.5. Accessories	18
	8.6. Instrument Settings	18
	8.7. Instrument Status	18
	8.8. Instrument Information	19
	8.9. Instrument Settings	19
~	o.to. Lump Settings	19
9.	USER ALCESS	20
	9.1. Adding d user	20
	9.3 Deleting a user	20
	9.4. Editing user access	20
10.	APPLICATIONS	21
	10.1. Single Wavelength	21
	10.2. Concentration via factor	22
	10.3. Wavescan	23
	10.4. KINELICS	24 25
	10.4.2 Parallel kinetics measurements	23
	10.5 Trace Manager – Overlauing & manipulating	21
	wavescan and kinetics files	29
	10.6. Standard Curve	31
	10.7. Equation Editor	33
11.	LIFE SCIENCE APPLICATIONS	39
	11.1. Nucleic Acid Applications	40
	11.1.1. DNA, RNA & Oligo	40
	11.1.2. CyDye™ DNA	42
	11.1.3. T <sub>m</sub> Calculation	43
	11.2. Protein Applications	46

11.2.1. BCA, Bradford, Lowry & Biuret Protein	
Assays	46
11.2.2. Determination of Protein Concentration	16
11.2.3. Determination of Protein Concentration	40
using direct UV methods	49
11.2.4. Protein UV	50
	51
12. SAVING AND PRINTING 12.1 Saving Sample Data	52 52
12.1.1. Internal	52
12.1.2. USB	52
12.1.3. USB csv	52
12.2. Automatic Saving	53
12.4. Exporting Data	53
13. SAMPLE MANAGER	54
13.1. Deleting data from the internal memory	54
13.2. Accessing Sample Manager from the main	
SCREEN 13.3 Accessing Sample Manager from within an	55
application	55
13.4. Recalled files	56
14. SAVING METHODS	57
14.1. Methods saved to the internal memory	57
14.2. Methods folder	57
14.5. Renaming methods tolder 14.4. Locking saved methods	58
14.5. Deleting saved methods	58
14.6. Backing up method folders to USB	59
14.7. Favourites folder	59
	59
15.1 Internal printer	60 60
15.2. Print via computer (PVC)	60
15.3. Bluetooth	60
15.4. Automatic printing	60
15.5. Manual printing	60
16.1 Operation	61 61
16.2. Saving Method & Results	62
16.3. Incorrect Fitting	62
16.4. Using the Cell Charger to Create a	~~
Standard Curve	62
17.1 Softing Sinner defaults	63
17.2. Calibration	64
17.3. Wash Mode	64
17.4. Cleaning/Maintenance/Troubleshooting	65
18. TECHNICAL SPECIFICATIONS	66
19. CHANGING TUNGSTEN AND DEUTERIUM LAMPS	67
20. TABLE OF ICONS	69
21. GLOSSARY OF BOXES	73

# Accessories Appendix

۱.	INSTALLATION	3
2.	<ul> <li>PASSIVE ACCESSORIES</li> <li>2.1. Single Cell Holder</li> <li>2.2. Variable Pathlength Cell Holder</li> <li>2.3. Water Thermostatted Cell Holder</li> <li>2.4. Film Holder</li> <li>2.5. Test Tube Holder</li> <li>2.6.1. General Information <ul> <li>2.6.1.1. General Warnings</li> <li>2.6.1.2. Quality, Reliability and Safety</li> <li>2.6.1.3. Safety Markings</li> </ul> </li> <li>2.6.2. Instrument Description <ul> <li>2.6.2.1. Controller front and rear panel</li> <li>2.6.2.2. Specifications</li> <li>2.6.2.2.1. Keypad/Display Unit</li> <li>2.6.2.2.3. Displayed Values</li> <li>2.6.2.4. Operating Specifications</li> <li>2.6.2.5. External Control Possibilities</li> <li>2.6.2.6. Dimensions</li> <li>2.6.2.7. Electrical data</li> <li>2.6.2.8. Description of Symbols</li> <li>2.6.2.9. Environmental conditions</li> <li>2.6.3.1. Checking the serial number</li> <li>4.6.3.2. Checking mains supply and fuses</li> <li>2.6.3.2.1. Changing the Mains Voltage Setting and Fuses</li> <li>2.6.3.3.1. Start up</li> <li>2.6.3.3.1. Start up</li> <li>2.6.3.3.2. External Connections</li> <li>2.6.4. Operating Instructions</li> <li>2.6.5.1. Cleaning the instrument</li> <li>2.6.5.2. Water level alarm</li> <li>2.6.6.1. Water level Alarm</li> <li>2.6.1.1. Water level Alarm</li> <li>2.6.1.2. Other system Alarms</li> </ul> </li> </ul>	4 5 5 5 5 5 6 6 6 6 7 7 7 7 7 7 7 7 7 7 7
3.	ACTIVE ACCESSORIES 3.1. 8-Position Cell Changer 3.2. 5-Position Thermostatted Cell Changer 3.3. Sipper & Flowcell 3.3.1. Sipper Troubleshooting 3.3.1.1. Sample not being sipped through Flowcell	11 11 12 12 13
	<ul><li>3.3.1.2. Sample being pumped right through Flowcell</li><li>3.3.1.3. Bubbles in Flowcell</li></ul>	14 14

3.3.1.4.	Inconsistent readings	14
3.3.1.5.	Wrong volume being sipped	14

15

4. INSTALLING THE INTERNAL PRINTER

# 1. HEALTH & SAFETY

### 1.1 Safety

There a number of warning labels and symbols on your instrument. These are there to inform you where a potential danger exists or particular caution is required. Before commencing installation, please take time to familiarize yourself with these symbols and their meaning.

This instrument is intended for use by individuals trained in and familiar with the use of spectrophotometers and their associated hazards. In the event of a malfunction or hazard occurring, the user responsible shall disconnect the unit from power and isolate for decontamination and /or repair.

This instrument is subject to the following hazards:



High voltages exist inside the unit. Repair and maintenance should only be carried out by individuals trained specifically to work on these instruments.



The UV source contained within the unit generates a light beam that traverses the sample chamber and is accessible in the lamp chamber. Under normal use the lamp beam is confined within the instrument. The unit should not be operated with the sample chamber lid open or the lamp housing lid removed. Prolonged exposure to the beam may cause permanent eye damage.



There are no bio-hazardous materials within the unit; however this unit could be used with bio-hazardous samples. Before using the instrument the customer should have in place decontamination procedures designed to protect laboratory workers from occupationally acquired infections. The sample chamber cell holders are removable and may be decontaminated using the appropriate disinfectant for the bio-hazard in question, rinsed with distilled water and then allowed to dry. The sample chamber and exterior may be wiped with a suitable disinfectant cleaning wipe.

- Decontamination. Equipment returned for repair should include an appropriate decontamination certificate.
- It is the responsibility of the customer to ensure that the user of the equipment is provided with a safe working environment.
- Any chemicals used in Analyses should be used, stored and disposed of in accordance with manufacturer's guidelines and local safety regulations.
- Toxic Fumes. Efficient laboratory ventilation must be provided when working with volatile solvents or toxic substances.
- Waste disposal. Disposal of some solvents and chemicals may be classed as hazardous waste and must be dealt with in accordance with local regulatory practice.
- Personal protective equipment. This is not required to operate the unit but the samples measured may require PPE. A local risk assessment should be carried out.



This equipment may be connected and controlled from a PC. To preserve the integrity of the measuring equipment it is essential that the attached PC itself conforms to basic safety and EMC standards and is set up in accordance with the manufacturers' instructions. If in doubt consult the information that came with your PC. In common with all computer operation the following safety precautions are advised.

- To reduce the chance of eye strain, set up the PC display with the correct viewing position, free from glare and with appropriate brightness and contrast settings.
- To reduce the chance of cross contamination from biological samples, use appropriate personnel protection measures and disinfectant wipes on keyboard and mouse.



Care must be taken when handling all heated accessories.

# 1.2. Unpacking & Installation

- These instruments weigh approximately 18.5 kg. Follow your local regulations for safe handling and lifting of this equipment.
- Inspect the instrument for any signs of damage caused during transit. If any damage is discovered. Do not use the instrument and report the problem to your supplier.
- The instrument must be placed on a stable, level bench or table capable of taking its weight with sufficient space around the instrument for ventilation to circulate freely.
- Ensure your proposed installation site conforms to the environmental conditions for safe operation.
  - Indoor use.
  - 5 to 40°C.
  - Maximum relative humidity 90% up to 31°C decreasing linearly to 50% at 40°C.
- Extremes of temperature may require re-calibration of the unit for optimum performance.
- If the instrument has been stored in a cold environment then it should be allowed to come to thermal equilibrium for 2 to 3 hours before operation so that start up calibration is not compromised by condensation.
- The equipment must be connected to the local supply outlet using the provided power cables. It can be operated from 90 to 264 V~ 50 or 60 Hz.
- Replace power inlet fuses only with the same type and rating as follows:

- For Deuterium/Tungsten units T 1.6 A H 250 V AC (Anti-Surge, High breaking capacity).
- For Xenon units T 1.6 A H 250 V AC (Anti-Surge, High breaking capacity).
- Power rating is:
  - 100 VA for Xenon units.
  - 150 VA for Deuterium/Tungsten units.
- The instrument should be positioned so that the power cable may be readily removed in the event of a hazard or malfunction occurring.
- Site the instrument in an atmosphere free from dust and corrosive fumes.
- Use the on/off switch on the left hand side of the instrument. The instrument will automatically perform some start up self diagnostic checks on switch on. Please wait for these to finish before attempting to use the equipment.
- 1.2.1. Instrument Connections





USB A plug for USB memory sticks (Ultrospec 7000, 8000 and 9000 stand alone units only)



Internal accessory connector

# 1.3. Equipment Operation

### 1.3.1. Controls and Indicators



### 1.3.2. Intended Use

The instrument is intended to be used by scientists and technicians who possess basic laboratory and technical skills and have the knowledge and understanding of the hazards involved, with the unit and the samples used, to operate it in a safe manner. e.g. DNA/RNA/Oligo concentration and purity measurements as well as protein determinations

### 1.3.3. Instrument Preparation

- Switch on the unit and allow it to finish its start up calibration.
- Best performance is obtained if the instrument is allowed to warm up and stabilize for at least 30 minutes.
- If applicable connect the unit to a PC using a USB cable and refer to the online help and user manual.
- Select the appropriate application or method.
- Where relevant, set up the application parameters for the sample.
- For Deuterium/ Tungsten units a precision mode is available. To use this mode enter the desired application and switch to precision mode.
- Select cuvette cells to use. It is important to use cells of the correct type. Most samples are measured using a standard 10 mm path length cell. Special cells and accessories are available for larger or smaller path lengths and sample volumes. It is important to use cells of the correct type. Some cells absorb UV and are not suitable for UV sample measurement. Cells used for samples should be free from dust, residue or scratches.

- Before preparing samples and sample reference blanks, you must be familiar with hazards arising from handling the sample materials and where necessary observe local regulatory practice, personnel protection equipment and measures designed to ensure your safety.
- Prepare the sample blanks (references). A reference is typically the solution that the sample is dissolved in. For Dual beam units the reference may be placed in the reference path or a separate sample reference may be made by first placing the reference in the sample path and performing a reference measurement.
- Prepare the sample solutions. The sample solution would normally comprise the sample under test dissolved in the reference solution.
- When placing the cells in the equipment ensure the cell is orientated so that the light energy will pass through the cell.
- For further information on running applications and methods refer to Chapter "10. *APPLICATIONS*".

### 1.3.4. Post Run procedures

- Empty cuvette cells of sample and rinse with deionized water.
- Clean cuvettes periodically with commercially available cleaning solution or dilute detergent solution followed by a thorough rinse in deionized water.
- Note that some samples and solvents may be classified as hazardous or bio hazardous waste. The disposal of such substances must be carried out in accordance with local regulatory practice.

### 1.3.5. Performance Validation

Good laboratory practice recommends that the unit is periodically checked for optical performance.

- Switch on validation checks. When the unit is powered up it performs wavelength accuracy and lamp energy checks. Once complete the unit will beep and show a green on light.
- Periodically wavelength, stray light and absorbance should be checked to ensure the unit is performing to specification. Deterioration in performance may indicate that the instrument needs servicing or that a poor baseline has been saved. Performance validation can be carried using traceable reference materials.

## 1.4. User Maintenance

- Other than the Tungsten and Deuterium lamps in certain instruments there are no user serviceable parts in this equipment.
- To prevent contamination:
  - Cell holders and accessories should be removed and cleaned with commercially available cleaning solution or dilute detergent followed by a thorough rinse in deionized water. Allow to dry thoroughly before use.
  - Casework and the sample compartment may be wiped down with commercially available disinfectant wipes.
- The lamps used in the unit age over time and less energy will be available to pass through the sample. Where energy has fallen to around 50% of the installation energy you are advised to have the lamps changed. Deuterium and Tungsten lamps may be replaced by the user, see section 20 for details. Xenon lamp replacement can only be carried out by a qualified service engineer.



If the equipment is operated in a manner not specified, then the protection provided by the equipment may be impaired and the instrument warranty withdrawn.

### 1.4.1. Troubleshooting

PROBLEM	POSSIBLE CAUSE
Start up calibration fails	Check sample (and reference) beam is clear before switch on. Possible optical failure, contact service support.
Negative absorbance reading	Check that sample and reference cells have not been swapped Check sufficient height of sample solution (beam height is nominally 15 mm above the cell base for normal measurements and it is recommended that it be filled to 20 mm above the base).
Unexpected results	Check for bubbles in solution.
Start up calibration fails	Check sample (and reference) beam is clear before switch on. Possible optical failure, contact service support.
Negative absorbance reading	Check that sample and reference cells have not been swapped.
	Check sufficient height of sample solution (beam height is nominally 15 mm above the cell base for normal measurements and it is recommended that it be filled to 20 mm above the base).
Absorbance values higher than expected	Check for use of incorrect cell type.
	Check cells are free from finger prints.
	Check cells for contamination.
	Check cell orientation.
	Check reference used.
	Possible optical alignment problem, contact service support.
Absorbance value lower than expected	Check sufficient height of sample solution.
	Check sample and reference are not the same.
	Check sample compartment lid is properly closed.
	Possible stray light issue, contact service support.
Instrument will not reference	Check orientation of cuvette.
	Check for use of incorrect cell type.
Poor reproducibility with nucleic acid	Check for particles, try using background correction.
anaiysis	ן וויט וווכופטצפ ווו כטווכפווגוטנוטוו.

# 1.5. Customer Support Contacts

**Note:** If you experience any problems with your instrument, please refer to section *"1.4.1. Troubleshooting"*.

If you require further assistance, please contact customer support :

visit:	www.biochromspectros.	com
or email:	support@hbiosci.com	
or telephone:	(US) 1-800-272-2775	(UK) +44 1223 423723
	1-508-893-8999	+44 1223 427888

### 1.6. Service, Repair or Return

Good laboratory practice recommends that the unit be periodically serviced to ensure the unit is suitable for purpose. It is recommended that the instrument be serviced annually. You can arrange this through your distributor. Prior to Inspection, Servicing, Repair or Return of Medical and Laboratory Equipment the unit must be decontaminated.

A returns policy operates on this equipment. Before returning the equipment to the distributor or manufacturer:

Please contact support@hbiosci.com for the correct paperwork and instructions.

## 1.7. Disposal

### Decontamination



In use this product may have been in contact with bio-hazardous materials. Before disposal all accessories should be removed and thoroughly cleaned in disinfectant and then rinsed with distilled water. All outside surfaces and sample chamber walls must be wiped down with disinfectant wipes suitable for purpose.

### WEEE



A label with a crossed-out wheeled bin symbol indicates that the product is covered by the Waste Electrical and Electronic Equipment (WEEE) Directive and is not to be disposed of as unsorted municipal waste. Any products marked with this symbol must be collected separately and in accordance with local regulatory practice.

# 2. INTRODUCTION TO THE Ultrospec 7000/8000/9000 SPECTROPHOTOMETER

The Biochrom Ultrospec 7000/8000/9000 are a range of standalone, easy to use, dual beam UV-visible spectrophotometers with high resolution color touch screens. All spectrophotometers offer a comprehensive range of spectrophotometric and life science applications.

A spectrophotometer is an optical arrangement that is designed to pass light energy through a sample. The reduction of transmittance of the light energy and corresponding absorbance peaks, caused by the sample under test, can be used in a qualitative way to determine sample composition or in a quantitative way, by comparing with known concentration standards, to determine the concentration of a sample. In other types of studies the tracking of absorbance over time can be useful to study chemical reactions and biological processes.

The Ultrospec 7000/8000/9000 range of spectrophotometers are designed to produce light energy from the far ultraviolet through visible light in the range 190 to 1100 nm. Many materials, and in particular solutions of materials will absorb light energy within this region. This makes the Ultrospec 7000/8000/9000 spectrophotometers applicable to a wide range of market sector needs including applications in Life sciences, clinical, pharmaceutical, cosmetics, food & drink, agricultural, industrial, environmental, toxicology, water treatment and teaching. There are a large number of published methods and assays available for these.

The icons described throughout this manual use the names quoted in Chapter "20. TABLE OF ICONS", if you are unsure of any of their functions please refer to this section.

For detailed descriptions of the functions of parameter boxes please refer to Chapter "21. GLOSSARY OF BOXES".

**Note:** Throughout this manual all screen shots are shown on a white background, this is purely for illustrative purposes.

# 3. USE WITH Datrys PC SOFTWARE

When connected to a PC the Ultrospec 7000/8000/9000 spectrophotometers can be controlled using the Datrys PC software package. Operation using Datrys PC software is described in the Datrys user manual or Datrys help file.

# 4. FREQUENTLY USED ICONS

Button	Name	Function		
٠	Forward arrow	Advances to the next screen in a sequence		
4	Back arrow	Returns to the previous screen in a sequence		
$\checkmark$	Tick	Confirms selection/entry. Saves and exits		
X	Cross/exit	Exit without saving		
Buttons	On The Sample	e Measurement Screen		
Ø	Take reference	Performs a reference measurement		
7	Take measurement	Performs a sample measurement		
<b>1</b>	Options arrow	Opens the options menu on the sample measurement screen		
Buttons On The Options Menu				
	Method parameters	Takes the user from the sample measurement screen to the first method parameter screen		
	Save data	Allows the user to manually save sample data to a specified location.		
0	Save method	Allows the user to save the current method parameters to the internal memory or a USB stick		
÷	Print	Prints the sample data from the specified printer		
	Auto print	Toggles auto print on and off		
	Sample Manager	Accesses Sample Manager		
4	Trace Manager	Accesses Trace Manager (wavescan and kinetics only).		

# 5. PERFORMING A MEASUREMENT

# 5.1. Ultrospec 7000/8000/9000

The Ultrospec 7000/8000/9000 are dual beam UV-Visible spectrophotometers that contain cell holders for both samples (situated at the front of the cell chamber) and references (situated at the back of the cell chamber). As dual beam spectrophotometers continually reference against the reference cell holder it is possible to perform measurements in one of three ways.

### To reference against air:

- **1.** Insert a cuvette containing the sample solution in the sample cell holder, ensure the reference cell holder is empty and close the sample chamber lid.
- 2. Press take measurement.
- **3.** Repeat until all sample data has been collected. See the section Saving and Printing for post measurement options.

### To reference against a solvent:

- **1.** Insert a cuvette containing the sample solution in the sample cell holder, insert a cuvette containing the solvent in the reference cell holder and close the sample chamber lid.
- 2. Press take measurement.
- **3.** Repeat until all sample data has been collected. See the section Saving and Printing for post measurement options.

Note: This methodology is used throughout the user manual.

### To correct for both solvent and cell effects (simulate matched cuvettes):

- **1.** Insert cuvettes containing the solvent in both the sample and reference cell holders and close the sample chamber lid.
- 2. Press take reference.
- **3.** Remove the cuvette containing solvent from the sample cell holder, empty out the solvent and replace with the sample solution.
- **4.** Insert the cuvette containing the sample solution in the sample cell holder and close the sample chamber lid.
- 5. Press take measurement.
- **6.** Repeat steps 4 and 5 until all sample data has been collected. See the section Saving and Printing for post measurement options.

Note: This methodology ensures the most accurate measurements.

# 6. LAMP MODE

When creating a method using an Ultrospec 7000/8000 or 9000 spectrophotometer the user will have the option of setting the Lamp Mode to Precision or Pulse. These modes are described below.

### Precision

Precision Mode is used for applications requiring the most accurate measurements. Both the Tungsten and Deuterium lamps will remain permanently switched on.

### Pulse

Pulse mode is used to help preserve lamp life and reduce running costs as the Tungsten and Deuterium lamps will be switched off after 15 minutes of inactivity.

**Note:** As the Ultrospec 7000 uses a Xenon lamp, measurements can only be made in pulse mode.

# 7. TYPES OF BOXES

Biochrom Ultrospec instruments use different kinds of boxes for parameter selection and entry. These include:

### Alphanumeric Text Entry

The alphanumeric text entry box allows the user to enter letters, numbers and symbols by pressing the abc, 123 and  $\lambda \mu$ ! buttons, respectively. It is possible to toggle between upper and lower case letters and through a list of symbols by pressing the abc and  $\lambda \mu!$  buttons twice.

Note: The layout of the screen is dependent on the text entry mode set under User Interface in Settings.



### Numeric Entry

The numeric entry box allows the user to include numbers in the method parameters. Depending on the numeric box selected it may be possible to add both positive and negative numbers.

Combo Box

Where there are greater than two options, the user will be presented with a combo box listing all of these. If there are greater than 8 options the user can scroll through these using either the page up and page down arrows or the scroll bar.

Note: If a box only contains two options i.e. On or Off, pressing this box will toggle between the options and not produce a combo box.





# 8. SETTINGS

Settings are accessed via the Settings button on the main screen (see below).



**Note:** When User Access Is turned on the Settings option will only appear on the main screen for users with Administrator or Supervisor privileges.



## 8.1. Date & Time

The Ultrospec 7000/8000/9000 will arrive with the US EST and date set. This can be changed by pressing the Date and Time button.

After the desired date and time have been entered, select the tick to save and exit or the cross to exit without saving.



## 8.2. Regional

The Ultrospec 7000/8000/9000 will arrive with the language set to English. This can be changed by pressing the Language button; the options are English, French, German, Spanish, Chinese and Japanese.

To save any alterations press the tick, to exit without saving press the cross.

Region	al
Language	
English	
Number Format	
999.9	
×	5

5061-046 Rev1

# 8.3. Data Output

This is the default saving and printing settings that will be used in all application method parameters.

### 8.4. User Interface

Allows the user to set the desired brightness level of the screen, the alphanumeric text entry mode and the duration after which the screensaver will be displayed (if required).

### 8.5. Accessories

Displays accessories that are fitted to the instrument and allows the user to set the desired accessory parameters. For specific details see the Accessories section.

# 8.6. Instrument Settings

The following options are included under instrument settings:

### 8.7. Instrument Status

This displays the current status of the instrument.





l

Peltier



Accessories

Cell Changer

1

Pathlength

Sipper



### 8.8. Instrument Information

Displays the serial number, user-interface (UI) version, build and release dates and instrument control (IC) version.

### 8.9. Instrument Settings

Instrument Settings allows the user to:

- **1.** Collect a new, temporary baseline. This will be stored until the instrument is switched off.
- **2.** Save the temporary baseline. This will become the permanent baseline and be stored until overwritten.
- **3.** Restore the original baseline. If measurements show the temporary baseline to be poor quality the permanent baseline can be restored.

**Note:** Using the Ultrospec 9000 the bandwidth at which the baseline is measured is selected by pressing the Bandwidth box

### 8.10. Lamp Settings

Displays the current lamp status and allows the user to reset lamp life (Tungsten and Deuterium lamps only).

To exit at any time press the exit button located in the bottom right hand corner.





Lamp Settings				
	Xen	on		
	Lamp State:	On		
	Installed Date: Installed Time:	1//1/2010 00:00:00		
	Original Energy: Current Energy:	20.7 20.7		
	Turned On: On Time: Total On Time:	108 9 9		
i	۹ 💽		×	

# 9. USER ACCESS

The Ultrospec 7000/8000/9000 have the option to assign users different access rights. These are set via the user access button.

Note: User access is only available to users who have administrator privileges.

Username	Password	Group
Administrator	1234	Admin
Alan	1234	Limited
Jo	1234	Superviso
Frank	1234	Limited
Helen	9999	Superviso
Steve	1234	Superviso
Katie	1234	Limited
Tarquin	1234	Limited

# 9.1. Adding a user

To add a new user to the instrument, press the 'Add user' button. The Ultrospec 7000/8000/9000 can store up to 16 individual users. Each user is given a user name (using alphanumeric entry), a 4-digit password and assigned to one of three user groups depending on the access level they require. The table below outlines the features each user group can access.

Password 1000 Group Base		
Password 1000 Group Base	ser Name	
1000 Group Base	Password	
Group Base	1000	
Base	Group	
	Base	

User Group	Run Applications & Saved Methods	Save Sample Data	Delete Sample Data from the instrument's	Save Methods	Access Settings Menu	Access User
			memory			Settings
Limited		$\checkmark$	×	×	×	×
Supervisor	1	<ul> <li>Image: A set of the set of the</li></ul>	$\checkmark$	<b>\</b>	×	×
Admin	1	<ul> <li>Image: A start of the start of</li></ul>	1			

### 9.2. Editing a user

To edit a user's details, highlight the desired user and press the 'Edit user' icon. This allows the username, password and user group to be edited/updated as above.

# 9.3. Deleting a user

To delete a user from the instrument, highlight the desired user and press the 'Add user' icon. Any methods or data created by this user will not be deleted. Note: It is not possible to delete the default administrator account.

# 9.4. Editing User Access

To disable user logins and user access, highlight the default administrator account and press the 'Edit User' icon to display the screen right.

**Note:** With Show Login set to No the instrument will not display the 'Switch User' button on the main screen and the instrument will always be in Administrator mode.



# **10. APPLICATIONS**



Single Wavelength	Absorbance, % transmission or concentration measurements at a single, specified wavelength.
Wavescan	Wavelength scan between two, user defined wavelengths in the range 190 to 1100 nm. The Ultrospec 7000/8000/9000 allows data overlay, post-scan data manipulation and user configurable peak and valley functions.
Kinetics	Serial and parallel measurements of absorbance versus time to determine rate or end points. The Ultrospec 7000/8000/9000 allows data overlay, post-scan data manipulation and user defined sectors.
Standard Curve	Concentration measurement at a single wavelength determined by the generation of a calibration curve of known standards.
Equation Editor	Allows users to create their own unique methods that include calculations and thresholds.

# 10.1. Single Wavelength

The Single Wavelength application performs simple absorbance (A) and % transmission (%T) measurements on samples, measuring the amount of light that has passed through a sample relative to a reference (this can be air).

### Measurement Parameters

Set Mode to Absorbance or %T. Wavelength, Bandwidth (Ultrospec 9000 only), Integration Time and Lamp Mode can be set as required. The Sample Seed entered under Sample will be the filename used for any data file saved automatically. You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.



Single Wavelength

 Auto-Print
 Auto Save

 On
 Off

 Print to...
 Save to...

 Internal Printer
 Internal

Set the outputs required in your method. For more information see the section Saving and Printing.

### **Taking A Measurement**

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.



### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

### 10.2. Concentration via factor

This mode within the Single Wavelength application makes simple concentration measurements on samples. Concentration is obtained by multiplying the measured absorbance at a specific wavelength by a factor. The factor may be known in advance or may be calculated by the instrument by measuring a standard of known concentration. Examples of concentration measurements include cholesterol, glucose and urea.

### **Measurement Parameters**

From the main screen of the Ultrospec 7000/8000/9000 select Applications followed by Single Wavelength to display the screen below.

Set Mode to Concentration. Wavelength, Bandwidth (Ultrospec 9000 only), Integration Time and Lamp Mode can be set as required.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.



Single Wavelength - Data Parameters
Factor Method
Predefined
Factor
324.1
Units
mg 100ml



Set Factor Method as required, enter Factor using numeric entry and Units using alphanumeric entry.

Set the outputs required in your method. For more information see the section Saving and Printing.

### **Taking A Measurement**

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

**Note:** For assays where there is no established concentration factor, calibration should be carried out using prepared standards (see Standard Curve for details of how to perform this).

### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

### 10.3. Wavescan

A measurement of the absorbance or %transmission of a sample over a specified wavelength range is one of the most useful physical characteristics of a compound, both as means of identification (qualitative analysis) and of estimation (quantitative analysis). The observed features arise due to the various electronic transitions that are possible within a molecule. The Ultrospec 7000/8000/9000 offer a range of post-scan data manipulation options including: 1st order derivative, enabling identification of multiple, unresolved peaks; 2nd order derivative, enabling identification of peak shoulders (inflections); 4th order derivative, which identifies both multiple peaks and inflections at the same time; Smoothing, utilizes the Savitzky-Golay algorithm to smooth data and increase the signal to noise ratio; Enhanced, which enhances features, sharpening peaks and valleys.

### Measurement Parameters.

Use Max Wavelength and Min Wavelength to set the required wavelength range (the Ultrospec 7000/8000/9000 scan from high to low wavelengths). Set Step, Mode and Scan Speed as required. If Scan Speed is set to Integration the user will also be required to set the Integration Time.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

Set Lamp Mode, Bandwidth (Ultrospec 9000 only) and Sample Overlays as required. The Sample Seed entered under Sample will be the filename of any automatically saved file.

**Note:** With Sample Overlays set to  $\geq 2$  the data all wavelength scans will be automatically saved to the instrument's internal memory and will be displayed in Trace Manager.







# 10.4. Kinetics

**Taking A Measurement** 

from the sample measurement.

Kinetics measurements made using a UV/visible spectrophotometer measure the change in absorbance at a single, fixed wavelength over a specified period. This can be used to provide useful information when an appropriate factor, defined in a reagent kit protocol, is applied. Reagent test kits are routinely used for the enzymatic determination of compounds in food, beverage and clinical laboratories.

UV/visible spectrophotometric kinetic assays are considered one of the most convenient measurements for enzymatic assays since they allow the rate of the reaction to be measured continuously.

This measurement parameters screen allows the user to set the following parameters:

Feature Detection: Determines the number of peaks or valleys that will be automatically detected. Options are Coarse, Sensitive or Custom.

Feature Type: The feature types that will be detected by the software. Options are Peaks or Valleys.

Feature Sort: Determines the how the features will be displayed in the data table. Options are Wavelength or Magnitude.

Draw Peaks: Switches the display of peak cursors on and off. The cursors show vertical dashed lines displaying the measured peak height and horizontal dashed lines showing the peak width.

Custom Peak Height: Only displayed if Feature Detection is set to Custom. This is the minimum height the peak has to be above the higher of the two adjacent minima for the peak to be detected.

Custom Peak Width: Only displayed if Feature Detection is set to Custom. This is the minimum width of the peak as determined by the difference in wavelength between the higher of the two adjacent minima and the opposing intersection of that higher minimum level and the peak profile.

Set the outputs required in your method. For more information see the section Saving and Printing.

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000)

insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take

screen will display a table below the scan. This table will display the Feature Type

selected in the method parameters. To manually add a peak or valley to the table,

position the cursor over the desired feature by either touching the feature or using

the left and right cursors and press an empty cell in the table.

#### Wavescan Feature Detection Draw Peaks Custom Off Feature Type Custom Peak Height Peaks 0.010 A Feature Sort Custom Peak Width Wavelength 50.00 nm





### 10.4.1. Serial Kinetics Measurements

Serial kinetics is the measurement of the absorbance of a single sample over a specified duration at a specified interval. As these instruments are capable of taking up to 1 reading per second serial kinetics measurements can be used for rapid rate reactions.

**Note:** Without a cell changer connected, the instrument is only capable of performing serial kinetics measurements.

### **Measurement Parameters**

*Wavelength, Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you require. The *Sample Seed* entered under Sample will be the filename of any automatically saved file. No. of Samples is described below.

**Note:** *Measurement Mode* only appears if a cell changer is connected. If this option is not displayed the instrument will perform a serial kinetics measurement.

**Note:** When performing a serial kinetics measurement with >1 sample the measurement proceeds as follows: Sample 1 will be measured for the full duration at the specified interval, after this is complete sample 2 will be measured for the full duration at the specified interval. The measurement will continue in this manner until all samples have been recorded. When measuring >1 sample, all data will be overlaid at the end of the measurement and automatically saved to the instrument's internal memory.

Set the *Delay* (time before first measurement), *Duration* (total measurement time, up to 180 minutes), *Interval* (duration between readings, from 1 second to duration) and *Integration Time* that you require.

**Note:** The integration time is determined by the interval (i.e. the maximum integration time is half the interval time).

*Mode* has options for Delta A, Final A and Slope and is the value that will be multiplied by the *Factor* to give the *Result* on the sample measurement screen. *Units* are entered using alphanumeric text entry and will appear on any printed or exported data. Y min and Y max are what is displayed during the measurement, the y axis auto-scales upon completion.







Set the outputs required in your method. For more information see the section Saving and Printing.

# Kinetics Print to... Auto Save Internal Printer On Auto-Print Save to... On Internal









### Taking a Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

Note: Measurements will commence after the specified *Delay* period (if applicable).

A measurement can be stopped at any time by pressing the Stop button at the bottom of the screen. All data collected to this point will be displayed on screen and can be saved.

The data displayed in the table below the scan refers to the full measurement range. To obtain data for a specific section it is necessary to add sections, this is done as follows: Set the cursor to the desired start position by either pressing on the scan or using the cursors, select t0 from the options menu, set the cursor to the desired end position and select t1 from the options menu. This can be repeated to add up to 4 discrete sections.

**Note:** Sections must be added in numerical order i.e. t1 must be added after t0, t2 after t1 etc.

With sections defined, the data displayed in the table below the scan is determined by the position of the cursor. Only when the cursor is positioned in the section of interest will this data be displayed.

A line of best fit can be added to any section by positioning the cursor in the desired section and selecting the add line of best fit button from the options menu.

**Note:** Details of how to perform overlays, data manipulation and selecting saved files can be found in the Trace Manager section.

### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

### 10.4.2. Parallel Kinetics Measurements

Parallel kinetics is the measurement of the absorbance of >1 sample (up to 8) at a set interval over a specified duration. E.g. for 4 samples a measurement would proceed as follows: The instrument will measure samples 1 to 4, pause for the appropriate interval, measure samples 1 to 4, pause and continue in this manner until the full duration has been reached.

**Note:** As parallel kinetics requires a cell changer the option of performing a parallel kinetics measurement is only available when a cell changer is connected.

### **Measurement Parameters**

*Wavelength, Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you require. The *Sample Seed* entered under *Sample* will be the filename of any saved file. No. of *Samples* is described above. *Measurement Mode* should be set to Parallel

Note: For parallel kinetics measurements the minimum number of samples is 2.

*Delay* and *Duration* (up to 180 minutes) can be set as required. The options available under *Interval* are determined by the number of samples set on the previous page.

Mode has options for Delta A, Final A and Slope and is the value that will be multiplied by the *Factor* to give the *Result* on the sample measurement screen. *Units* are entered using alphanumeric text entry and will appear on any printed or exported data. Y min and Y max are what is displayed during the measurement, the y axis auto-scales upon completion.

Set the outputs required in your method. For more information see the section Saving and Printing.









As parallel kinetics requires a cell changer, you will be given the option of viewing the prompt for sample position before collecting data. All other cell changer parameters are determined by the method parameters. This prompt is shown below.











### Taking a Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000), insert a cuvette containing the reference solution in the reference cell holder and a cuvettes containing samples in positions 1 through 8 and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

A measurement can be stopped at any time by pressing the Stop button at the bottom of the screen. All data collected to this point will be displayed on screen and can be saved.

The data displayed in the table below the scan refers to the full measurement range of the selected trace (see below for selecting a trace). To obtain data for a specific section it is necessary to add sections, this done as follows: Set the cursor to the desired start position by either pressing on the scan or using the left and right cursors, select t0 from the options menu, set the cursor to the desired end position and select t1 from the options menu. It is possible to add up to 4 discrete sections.

**Note:** Sections must be added in numerical order i.e. t1 must be added after t0, t2 after t1 etc.

With sections defined, the data displayed in the table below the scan is determined by the position of the cursor. Only when the cursor is positioned in the section of interest will this data be displayed. This refers to the selected scan only and is indicated by the color of the column headers in the table.

A line of best fit can be added to any section by positioning the cursor in the desired section and selecting the add line of best fit button from the options menu.

To allow post scan data manipulation all samples measured are automatically included in Trace Manager (see below for full details of Trace Manager).

**Note:** The data displayed in the table below the scan refers to the <u>Selected</u> sample only.





### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

### 10.5. Trace Manager – Overlaying & Manipulating Wavescan & Kinetics Files

Trace Manager is the application used by the Ultrospec 7000/8000/9000 to overlay and manipulate wavescan and kinetics files. Samples are loaded into Trace Manager as described below:

### Using Sample Manager

Wavescan and kinetics files can be loaded directly into Trace Manager by selecting the required files from Sample Manager on the main screen. This procedure is outlined below:

Highlight the required files by pressing on the appropriate row and load these into Trace Manager by pressing the Sample Manager icon (bottom right corner). If the files you require are not displayed on the screen you can scroll through the pages using the up and down arrows. Pressing on the Sample ID and Date column headers will sort the files alphabetically and chronologically, respectively.

**Note:** If a USB memory stick is inserted, it is possible to toggle between files stored on the internal and USB memories using the icon in the left hand corner. The icon in the top right corner of the screen will display the memory that is currently in use.

All recalled files will be displayed on the Trace Manager screen (up to 8). The color of the *Slot* number will be the color of the trace when it is displayed on screen. To toggle which files are displayed on the measurement screen press in the appropriate *Display* box (up to 8 files can be displayed). To select which files data will be displayed on the measurement screen press in the appropriate *Selected* box (only one file may be selected). Press the forward arrow to view the overlaid data.

Sample Id	Application	Date	4
Tank1	Wavescan	2010/07/09 08:33:22	Ĩ
Tank2	Wavescan	2010/07/09 08:33:24	
Tank3	Wavescan	2010/07/09 08:33:25	
Tank4	Wavescan	2010/07/09 08:33:27	
Tank5	Wavescan	2010/07/09 08:33:29	ĺ
Tank6	Wavescan	2010/07/09 08:33:31	
Tank7	Wavescan	2010/07/09 08:33:34	Ĭ.
Tank8	Wavescan	2010/07/09 08:33:36	

		Trace Manag	ger	
Slot	Source	Trace Output	Display	Selected
1	Tank1	Sample Data	1	1
2	Tank2	1st Derivative	1	×
3	Tank3	2nd Derivative	1	×
4	Tank4	4th Derivative	1	×
5	Tank5	Smoothed	1	×
6	Tank6	Enhanced	1	×
7	Tank7	Sample Data	1	×
8	Tank8	Sample Data	1	×
			â	× •

5061-046 Rev1

### From Within An Application

To overlay saved files with a live trace displayed, Trace Manager can be accessed from within the application. This procedure is required for Limited users as they do not have the ability to access Sample Manager on the main screen.

Trace Manager can be accessed from the Wavescan and Kinetics measurement screens using the Trace Manager icon on the options menu.

When accessed with no overlays displayed, Trace Manager will be empty. Files are added to this screen by selecting the Sample Manager icon in the left hand corner of the screen and loading saved files as described below.

Highlight the required files by pressing on the appropriate row and load these into Trace Manager by pressing the Sample Manager icon. If the files you require are not displayed on the screen you can scroll through the pages using the up and down arrows. Pressing on the Sample ID and Date column headers will sort the files alphabetically and chronologically, respectively.

**Note:** If a USB memory stick is inserted, it is possible to toggle between files stored on the internal and USB memories using the icon in the left hand corner. The icon in the top right corner will display the memory that is currently in use.

All recalled files will be displayed on the Trace Manager screen (up to 8). The color of the *Slot* number will be the color of the trace when it is displayed on screen. To toggle which files are displayed on the measurement screen press in the appropriate *Display* box (up to 8 files can be displayed). To select which files data will be displayed on the measurement screen press in the appropriate *Selected* box (only one file may be selected). Press the forward arrow to view the overlaid data.

### Post Scan Manipulation

Trace Manager allows the user to manipulate recalled wavescan and kinetics data using the procedure outlined below.



		Trace Manag	jer	
Slot	Source	Trace Output	Display	Selected
1				
2				
3				
4				
5				
6				
7				
8				
			â	× +

Sam	ple Manager -	page 1 of 4	
Sample Id	Application	Date	۵
Tank12	Wavescan	2010/02/09 08:32:35	
Tank11	Wavescan	2010/02/09 08:32:33	۵
Tank10	Wavescan	2010/02/09 08:32:32	۵
Tank9	Wavescan	2010/02/09 08:32:30	
Tank8	Wavescan	2010/02/09 08:32:29	۵
Tank7	Wavescan	2010/02/09 08:32:27	
Tank6	Wavescan	2010/02/09 08:32:26	
Tank5	Wavescan	2010/02/09 08:32:25	
• <del>~~</del> +	<b>.</b>	Х ≙ ◆	

		Trace Manag	ger	
Slot	Source	Trace Output	Display	Selected
1	Tank1	Sample Data	1	1
2	Tank2	1st Derivative	1	×
3	Tank3	2nd Derivative	1	×
4	Tank4	4th Derivative	1	×
5	Tank5	Smoothed	1	×
6	Tank6	Enhanced	1	×
7	Tank7	Sample Data	1	×
8	Tank8	Sample Data	1	×
			â	× +

With the required files loaded into Trace Manager press the appropriate *Trace Output* box to display the manipulation options.

		Trace Manag	jer	
Slot	Source	Trace Output	Display	Selected
1	Tank1	Sample Da	ita	1
2	Tank2	Sample Da	ta	×
3	Tank3	1st Derivat	ive	×
4	Tank4	2nd Derival	iue	×
5	Tank5			×
6		4th Derivat	ive -	
7		Smoothee	4	
8		Enhance	a –	
			â	× +

Wavescan post scan manipulations are:

Sample Data	Displays the raw wavescan data (this is the default option).
1st – 4th Derivative	Displays the derivative data to the desired order.
Smoothed	Uses the Savitzky-Golay algorithm to reduce noise and smooth the data.
Enhanced	Enhances features, sharpening peaks and valleys.

Kinetics post scan manipulations are:

Sample Data	Displays the raw kinetics data (this is the default option).
Low	Applies a low level of smoothing to the data.
Medium	Applies a medium level of smoothing to the data.
High	Applies a high level of smoothing to the data.

After the manipulations and display options have been set as required press the forward arrow in the right corner to display the recalled data on the sample measurement screen.

**Note:** For wavescan and serial kinetic measurements with overlays set to >1 and parallel kinetic measurements, data will be automatically saved into the instrument's internal memory and displayed in Trace Manager. To ensure the optimum performance of the instrument it is recommended that unwanted files are deleted from the internal memory at regular intervals (see Sample Manager – deleting files from the internal memory).



# 10.6. Standard Curve

The construction of a multi-point calibration curve from standards of known concentration to quantify unknown samples is a fundamental use of a spectrophotometer. The Ultrospec 7000/8000/9000 range of instruments has the advantage of being able to store calibration curves with a method. Each calibration curve can be created using up to 9 standards, with each standard measurement being made of up to 3 replicates.

### Creating a Standard Curve

Set *Wavelength, Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* as required. The Sample Seed entered under *Sample* will be the filename of any file saved automatically.



*Calibration can be set to Standards* (the user is required to prepare and measure standards) or *Manual* (the user inputs both the standard concentrations and standard absorbances). Set *Standards, Replicates, Curve Fit* and *Units* as required in your application.

Set the concentration values for each of the standards using the numeric entry box.

Set the outputs required in your method. For more information see the section

To create the standard curve when using replicates press the replicates button in the bottom right corner to take you to the screen shown overleaf. With Replicates off standards can be measured directly as described overleaf.

**Note:** Pressing the save method icon before any standards have been measured will save the method parameters only. Recalling a method containing method parameters only requires the user to construct a standard curve before measuring samples.

To create a standard curve using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the first standard/replicate in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the standard measurement.

Continue recording all standards/replicates until the standard curve has been completed. To repeat any standard measurements simply press the desired result, insert the correct standard and press take measurement.

**Note:** After all replicates have been taken for a particular standard pressing the replicates icon takes the user to the next standard that was specified in the method.

Saving and Printing.











To ignore any outlying standard measurements, press the tick in the appropriate row to toggle it to a cross. Any ignored measurement will be automatically removed from the standard curve. These can be reinstated by pressing the cross.

#### Standard Curve Standards 0.25 2.500 0.050 A × 0.20 1 3.000 0.100 A 0.15 x 3.500 0.150 A 1 0.10 0.200 A 4.000 1 4.500 0.250 A = 0.0492x, F У ø A

### **Taking A Measurement**

After the standard curve has been collected press the forward arrow to proceed to the sample measurement screen.

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

To view the standard curve whilst on the sample measurement screen, simply press the 'View Curve' icon that appears under the options menu.

**Note:** Saving the method using the Save Method icon that appears on the options menu will save both the method parameters and the standard curve. Recalling a method containing method parameters and standard curve allows the user to measure samples directly.

### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

### 10.7. Equation Editor

The Equation Editor application allows users to create their own unique methods that include calculations and thresholds. Examples of methods that can be created using Equation Editor include percentage strength calculations and olive oil and chlorophyll analyses.

### **Getting Started**

The use of Equation Editor is outlined below.

#### **Measurement Parameters**

The first screen of Equation Editor allows the user to set the measurement parameters that will be used for all subsequent measurements. *Bandwidth* (Ultrospec 9000 only), *Integration Time, Lamp Mode* and *Mode* are used as in all other applications. The use of *Prompt between*  $\lambda$  and Scan are described below.







speeds up analyses as only the required wavelengths are measured.

Scan on The instrument will perform a full wavelength scan over a range set under the sample measurement parameters screen. With scan on the user can set both the scan speed and the step size.

You can advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

Using the Sample Measurement table the user inputs all of the measurements that are required in the method i.e. the screen left shows a fixed wavelength measurement at 430.0 nm that is named 430. Data is inputted as described on next page. Before any data is inputted Name will read 'Not set' and will not appear in the Sample Data list on the Equation Builder.

Equation Editor - Samples				
Name	λ	Function	Limit	Del
430	430.0	Abs at λ		×
Not set	0.0			×
Not set	0.0			×
Not set	0.0			х
Not set	0.0			×
Not set	0.0			×
Not set	0.0			×
Not set	0.0			×
4	•		•	
•	-		7	

- Name Pressing this allows the user to input an alphanumeric name for the measurement. The name inputted here will be used as the Sample Data in the Equation Builder (inputted data only appears in the Equation Builder if the user has defined a name).
- Wavelength Pressing this produces the number entry box that allows the user to input the ( $\lambda$ ) wavelength for measurement.

Function Pressing this displays a combo box with the following options.

Abs / %T at  $\lambda$  – measurement will be at the wavelength inputted by the user only.

Valley closest to  $\lambda$  – the instrument automatically finds the valley closest to the inputted wavelength.

- +/- Used with Peak closest to  $\lambda$  and Valley closest to  $\lambda$  only. This is the range over which the instrument will scan for a peak or valley from the inputted wavelength.
- Del Pressing this deletes the row. If the data is used in an equation this will also be deleted.

Equation Editor has four options for sample naming conventions. These are shown below:



Default	The sample name consists of Sample and an incrementing number.
Auto Increment	The sample name is a combination of sample seed and an incrementing sample number. The user will be prompted to enter the sample seed for each new batch of samples.
Prompt for ID	The user is prompted to enter the sample name before running each sample.
Fixed List	The user will be prompted to enter the number of samples required. Sample names for each sample are then entered on the subsequent screens. The inputted sample names are saved for each method.

The Standard Specification screen is used to declare a list of all of the standard solutions which will be referenced when creating an equation.

E.g. the equation for percentage strength at 500 nm compares the absorbance of a sample to the absorbance of a control standard. This is where the control standard would be defined.

Standard names are entered by pressing on the desired row and entering the standard name using alphanumeric text entry. Standard measurements can be made for any measurements specified in the Samples table.

The Constant Factor Specification screen is used to declare any constants used in the equation. Data is inputted as described below. Before any data is inputted *Constant Name* will read 'Not set' and will not appear in the Constants list on the Equation Builder. Constant Name Pressing this allows the user to input an alphanumeric name for the constant

Value	Pressing this produces the number entry box that allows the user
	to input the value of the constant.

Units Pressing this allows the user to enter units for the constant using alphanumeric text entry. If this column is left blank, no units will be displayed on exported or printed data.

Del Pressing this deletes the row. If the constant is used in an equation this will also be deleted.

The Variable Factor Specification screen is used to declare any variables used in the equation. Data is inputted as described below. Before any data is inputted, *Variable Name* will read 'Not set' and will not appear in the Variables list on the Equation Builder.

		L
Variable Name	Pressing this allows the user to input an alphanumeric name for the variable factor.	
Default	Pressing this produces the number entry box that allows the user to input the default variable factor. Default values can be edited during a measurement.	
Jnits	Pressing this allows the user to enter units for the variable factor using alphanumeric text entry. If this column is left blank, no units will be displayed on exported or printed data.	
Change Each	Pressing this toggles between sample and batch. When set to sample the method will prompt for a variable to be entered before each sample measurement. When set to batch, the method will prompt for a variable to be entered at the start of each sample batch	

Equation Editor - Standard Specification	1
Standard Names	Del
Standard 1	х
Not set	х
Not set	х
Not set	×
Not set	×
Not set	х
Not set	×
Not set	×
• •	

Equation Editor - Constant Factor Specification				
	Constant Name	Value	Units	Del
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	Not set	1.000		×
	4		•	
	T			

Equation Editor - Variable Factor Specification				
Variable Name	Default	Units	Change On	Del
Not set	1.000		Sample	×
Not set	1.000		Sample	х
Not set	1.000		Sample	×
Not set	1.000		Sample	×
Not set	1.000		Sample	х
Not set	1.000		Sample	×
Not set	1.000		Sample	х
Not set	1.000		Sample	×
4			•	
•			-	

The Equation Viewer screen provides an overview of any equations created, as well as allowing the user to create new or edit existing equations. Data is inputted as described below. Before any data is inputted Name will read 'Not set' and will not appear on the results screen.

on Editor - Equation V	√iewer	
Equation	Units	Del
	*	х
		х
		×
		х
		×
		×
		х
		х
	•	
	n Editor - Equation Equation	Equation Viewer Equation Units   Equation

- Name Pressing this allows the user to input a unique, alphanumeric name for the equation. The names entered here will be displayed in the 'Equations' combo box on the Equation Builder screen as well the results screens. Therefore any equation can be easily identified when using it in other equations.
- Equation Pressing this takes the user to the 'Equation Builder' (see below). Any equation constructed in the Equation Builder will be displayed in this box.
- Units Pressing this allows the user to enter units for the result of the equation using alphanumeric text entry. If this column is left blank no units will be displayed alongside the result.
- Del Pressing this deletes the row and removes the equation.

The Equation Builder allows the user to create any equations required in the method. Data is inputted as described below. To allow data to be inserted or deleted the cursor can be moved left and right using the left and right arrows. Data can be deleted using the delete icon and thresholds added using the thresholds icon (see below).



Variables	Pressing on this displays a list that contains any variables added to the 'Variables' table by the user. Selecting the desired variable enters it into the equation.
Constants	Pressing on this displays a list that contains any constants added to the 'Constants' table by the user. Selecting the desired constant enters it into the equation.
Equations	Pressing on this displays a list that contains any equations that have been created in this method. Selecting the desired equation enters it into the equation.
Sample Data	Pressing on this displays a list that contains all of the readings specified by the user in the Sample Measurement table. Selecting the desired sample data enters it into the equation.
Symbols	Pressing on this displays a list containing mathematical symbols and the logic gates AND and OR. Selecting the desired symbol enters it into the equation.
Numbers	Pressing this produces the number entry box that allows the user to directly input numbers into the equation.
Note: All of th	a data above appears in the lists as it was inputted in the appropriate

**Note:** All of the data above appears in the lists as it was inputted in the appropriate table.

After the equation has been inputted the user has two options. If the result is to be viewed as a number, press the back arrow to return to the Equation Viewer screen. If the result is to be viewed with a user specified pass/fail limit press the 'Thresholds' button to set appropriate thresholds for the measurement.
The first thresholds screen allows the user to input how many thresholds are required for a result.





With Thresholds set to 3 the screen will display the table layout shown left. Setting Thresholds to 2 and 1 will reduce the number of values you can input to 2 and 1, respectively.

- Value Pressing this produces the number entry box that allows the user to directly input numbers for threshold values.
- Name Pressing this allows the user to enter the text that will be displayed for each result using alphanumeric entry.

Using the example above: A result of  $\geq 100$  will return the answer Position 1, a result of <100 and  $\geq 50$  will return the answer Position 2, a result of <50 and  $\geq 10$  will return the answer Position 3 and a result of <10 will return the answer Position 4.

**Note:** When using AND or OR logic in an equation the result will be returned as a binary (1 = true and 0 = false). These can be incorporated into the thresholds by setting the Thresholds to 1, Value to 1 and setting the appropriate names (as the name above value corresponds to  $\geq 1$  this is the response that will be displayed for true results).

After all of the thresholds have been set, pressing the forward arrow will return the user to the Equation Builder screen. The Equation Viewer screen can be accessed using the back arrow.

Continue as described above until all equations have been created. Once complete pressing the forward arrow will take the user to the output options screen below.

Set the outputs required in your method. For more information see the section Saving and Printing.

**Note:** Automatically saving sample data with sample naming set to 'Prompt for ID' or 'Fixed list' will save the data using the first inputted sample name as the filename.

Equatio	n Editor
Print to	Auto Save
Internal Printer	On
Auto-Print	Save to
On	USB
•	•

5061-046 Rev1

#### Performing A Measurement - No Standards

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

**Note:** Pressing on the equation name allows the user to view the equation and the individual measurement results.

## Using Standards

When performing a measurement using a method that includes standards, the first press of the take measurement button will produce a message box that prompts the user to insert a specific standard. After all standards have been measured, subsequent presses of the take measurement button will perform sample measurements.

#### Saving Methods And Exiting

As methods developed using Equation Editor may have taken time to input, the software will prompt the user to save the method before exiting. Pressing the cross will return to the user to the results screen, where they can save the method, pressing the tick will exit without saving. Details of method saving can be found in the Saving Methods section.

#### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

	I	Equatio	n Edita	r		
		Sar Sam	n <b>ple</b> ple1			
Equ	ation		Resu	ilt	Un	its
% (	ourity		98.	5	2	6
B	esult		Pass	5		
			ŧ	×	Ē	





# 11. LIFE SCIENCE APPLICATIONS

This contains two sub folders; Nucleic Acids and Protein. The contents of these sub folders are detailed below:



#### **Nucleic Acids**

inaciere neras	
DNA	Utilizes the absorbance measurements at 230, 260 & 280 nm with optional background correction to perform a concentration and purity check for DNA samples.
RNA	Utilizes the absorbance measurements at 230, 260 & 280 nm with optional background correction to perform a concentration and purity check for RNA samples.
Oligo	Utilizes the absorbance measurements at 230, 260 & 280 nm with optional background correction to perform a concentration and purity check for oligo samples.
CyDye DNA	Measures the labelling efficiency of fluorescently labelled DNA probes to ensure that there is sufficient amount of each probe to give satisfactory signals. The DNA yield is measured at 260 nm whilst the incorporation of the dyes is measured at the absorption maxima. This method is also useful for measuring the yields and brightness of fluorescently labelled in-situ hybridization probes.
T <sub>m</sub> Calc	The $T_m$ Calculation application calculates the theoretical melting point from the base sequence of a primer. It is done using nearest neighbour thermodynamic data for each base in the nucleotide chain in relation to its neighbour.
Protein	
BCA™	Quantitative determination of protein concentration utilising the absorbance measurement at 562 nm.
Bradford	Quantitative determination of protein concentration utilising the absorbance measurement at 595 nm.
Lowry	Quantitative determination of protein concentration utilising the absorbance measurement at 750 nm.
Biuret	Quantitative determination of protein concentration utilising the absorbance measurement at 546 nm.
Protein UV	Direct UV determination of protein concentration at 280 nm using the Christian Warburg calculation.
Protein A280	Direct UV determination of protein concentration using BSA, IgG, Lysozyme, Molar Extinction. Mass Extinction or E1% calculations.

## 11.1. Nucleic Acid Applications

## 11.1.1. DNA, RNA & Oligo

Nucleic acids can be quantified at 260 nm because it is well established that solutions of DNA and RNA in 10 mm pathlength cells with an optical density (absorbance) of 1.0 have concentrations of 50  $\mu$ g/ml and 40  $\mu$ g/ml, respectively. Oligonucleotides typically have a factor of 33  $\mu$ g/ml, although this does vary with base composition and can be calculated if the base sequence is known.

### Concentration = $A260 \times Factor$

Ultrospec 7000/8000/9000 use the default factors 50, 40 and 33 for DNA, RNA and oligonucleotides, respectively. Compensation for dilution and pathlength can also be entered.

## Nucleic Acid Purity Checks

- Nucleic acids extracted from cells are accompanied by proteins and extensive purification is required to separate these protein impurities. The ratio of A260/A280 gives an indication of a sample's purity, with pure DNA and RNA preparations typically having ratios of ≥1.8 and ≥2.0, respectively. Deviations from these values indicate the presence of impurities, but care must be taken when interpreting results.
- Concentration also affects both the A260 and A280 readings. If a solution is too dilute, the readings may be at the instrument's detection limit and results may vary as there is less distinction of the A260 peak and A280 slope from the background absorbance. For accurate measurements A260 should always be greater than 0.1.
- Elevated A230 values can also indicate the presence of impurities (230 nm is near the absorbance maximum of peptide bonds and also indicates buffer contamination since EDTA and other buffer salts absorb at this wavelength). When measuring RNA samples, the A260/A230 ratio should be >2.0. Ratios lower than 2.0 generally indicate contamination with guanidinium thiocyanate, a reagent commonly used in RNA purification and which absorbs over the 230-260 nm range. A wavelength scan of the nucleic acid is particularly useful for RNA samples.

## **Background Correction**

- To compensate for the effects of background absorbance caused by turbidity, high absorbance buffer solutions and the use of reduced aperture cells the Ultrospec 7000/8000/9000 have the option of background correction at a 320 nm.
- When used, A320 is subtracted from A260 and A280 prior to use so that: Concentration = (A260 - A320) × Factor Abs ratio = (A260 - A320) / (A280 - A320) Abs ratio = (A260 - A320) / (A230 - A320)
- The use of background correction can remove variability due to handling effects of low volume disposable cells.

### Spectral scan of nucleic acid

**Note:** An absorbance maximum near 260 nm and absorbance minimum near 230 nm, a flat peak near 260 nm and steep slope at 280 nm and very little absorbance at 320 nm.



#### **Measurement Parameters**

Set *Pathlength* and *Dilution Factor* to the required values. Check above to see if Background correction is required. To perform a scan over the range 220–330 nm for each measurement set *Scan* to *On*. Set Units to encompass the expected concentration of your samples, the default *Factor* will update automatically depending on the units selected i.e. for units of  $\mu$ g/ml the default factor will be 50.00. If the Factor required differs from the default value this can be edited using numeric entry.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

*Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you require. The *Sample Seed* entered under *Sample* will the filename of any saved file.

Set the outputs required in your method. For more information see the section Saving and Printing.







## Taking A Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

If *Background* is set to *On* the A320 result will be included in the left hand column and automatically subtracted from the displayed A230, A260, A280, A260/A230, A260/A280 results.

If *Scan* is set to *On* in method parameters a scan of the most recently run sample can be viewed by pressing the View Scan icon in the options menu.

**Note:** When saving sample data, scan files will not be saved. The *Wavescan* application should be used to save scans of nucleic acid samples.





#### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

## 11.1.2. CyDye DNA

The measurement of the labelling efficiency of fluorescently labelled DNA probes before 2-color micro-array hybridization ensures that there is sufficient amount of each probe to give satisfactory signals. The data also provides an opportunity to balance the relative intensities of each fluorescent dye by adjusting the concentration of each probe before hybridization. The DNA yield is measured at 260 nm whilst the incorporation of the dyes is measured at the absorption maxima. This method is also useful for measuring the yields and brightness of fluorescently labelled *in-situ* hybridization probes.

## **Measurement Parameters**

Number of Dyes, this can be set to 1 or 2. Dye 1 Name allows the user to select the dye used in the measurement, the Ultrospec 7000/8000/9000 have 19 dyes pre-programmed and the option for user entry using the Custom Dye option.  $\lambda$ *Max, Extinction Coefficient* and *Correction Factor* are only editable when using the Custom Dye option.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

With *Number of Dyes* set to 2, the next method parameter screen allows the user to specify the second dye used in the measurement.

If the measurement requires the calculation of DNA Concentration and/or DNA Quantity the relevant nucleic acid can be inputted. A custom factor can be entered by selecting Custom in *Nucleic Acids*.









This parameters screen allows the user to set the pathlength of the cuvette being used, whether background correction is required and at what wavelength, the dilution factor and volume of the sample in  $\mu$ l. All of these will be used in the calculations.

*Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you wish. The *Sample Seed* entered under Sample will the filename of any saved file.

Set the outputs required in your method. For more information see the section Saving and Printing.

# CyDye DNA Auto-Print Auto Save On Off Print to... Save to... Internal Printer Internal





### Taking a Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

If *Background Correction* is set to *On* the *Background Wavelength* set in the method will be included in the left hand column and automatically subtracted from all results.

To toggle between DNA & dye parameters press the toggle icon on the options menu. CyDye DNA measurements are wavelength scanning measurements, to view the scan for a particular measurement press the scan button in the options menu.

## Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

## 11.1.3. T<sub>m</sub> Calculation

The  $T_m$  Calculation application calculates the theoretical melting point from the base sequence of a primer. It is done using nearest neighbour thermodynamic data for each base in the nucleotide chain in relation to its neighbour (Breslauer *et al*, *Proc.Natl. Acad. Sci. USA*, **1986**, 83, 3746). The data obtained are useful in both the characterization of oligonucleotides and in calculating  $T_m$  for primers used in PCR experiments.

The ACGT/U sequence entered in the method parameters is used to calculate the theoretical  $T_m$ , the theoretical absorbance (Absorbance units/mmol) and the conversion factor (mg/ml). This is possible as the stability of a bent and twisted sequence of bases such as an oligonucleotide is dependent on the actual base sequence. These calculated thermodynamic interactions between adjacent base pairs have been shown to correlate well with experimental observations.

The  $\rm T_m$  Calculation application uses matrices of known, published thermodynamic values and extinction coefficients to calculate  $\rm T_m$  and the theoretical absorbance/factor of an entered base sequence.

 $T_m$  is calculated using the equation:-

 $T_{m} = \frac{\Delta H \times 100}{\Delta S + (1.987 \times \log_{e}(c/4 + 53.0822))} - 273.15 + \log [salt]$ 

where  $\Delta H$  and  $\Delta S$  are the enthalpy and entropy values, respectively summed from respective 2 × 4 × 4 nearest neighbour matrices

c is the Primer concentration of oligonucleotide (pmoles/ml) in the calculated  $T_m$  or the measured concentration in measured  $T_m$ . In the latter case, concentration is obtained from the equation:

c = Abs(260 nm) × Calculated factor × pathlength multiplier × 10 000

MW

Calculated factor and MW are defined below.

**[salt]** is the buffer molarity plus total molarity of salts in the hybridization solution (moles/l).

Weights for  $\Delta S$  are indexed by adjacent paired bases. A similar equation applies to weights for  $\Delta H$ , again indexed by adjacent bases.

**Note:** Bivalent salts may need normalising using a multiplying factor of 100 because of their greater binding power.

## **Theoretical Absorbance**

The Theoretical Absorbance is based on a calculation as follows:

For each adjacent pair of bases (nearest neighbours) an extinction coefficient weight is accumulated using a  $4 \times 4$  table (one for either DNA or RNA). This total weight is doubled and then for each internal base a counterweight is subtracted using another  $1 \times 4$  table. The end bases are excluded from the latter summation.

Total Extinction Coefficient E =  $\frac{\sum (2 \times aTable[base\_type][base(n)][base(n+1)])}{\sum (tTable[base\_type][base(n]])}$ 

## **Conversion Factor**

The Conversion Factor is given by =

 $\frac{\text{Molecular weight}_{\text{ABCDE}}}{\Sigma_{\text{EABCDE}}}$ 

where

 $E_{ABCDE} = [2 \times (EAB + EBC + ECD + EDE) - EB - EC - ED]$ 

• The molecular weight (MW) of a DNA oligonucleotide is calculated from:

 $\mathbf{MW} (g/mole) = [(dA \times 312.2) + (dC \times 288.2) + (dG \times 328.2) + (dT \times 303.2.]] + [(MWcounter-ion) \times (length of oligo in bases)] (for RNA oligonucleotide, (dT \times 303.2) is replaced by (dU \times 298.2))]$ 

The MW calculated using this equation must be adjusted for the contribution of the atoms at the 5' and 3' ends of the oligo.

For phosphorylated oligos:	Add $[17 + (2 \times MW \text{ of the counter-ion})]$
For non-phosphorylated oligos:	Subtract [61 + (MW of the counter-ion)]

The MW (g/mole) of the most common oligo counter ions are:

Na (so	dium)	23.0
K (potc	issium)	39.1
TEA (tri	ethylammonium) 102.2	
Other	Defaults to 1.0 (variable	0.1-999.9)

Calculated molecular weight: a weight is added for each base looked up from a table. The weight of the counter ion is added for every base from a small table for the known ions. If phosphorylated, then the system adds 17.0 plus two counter ions otherwise it subtracts 61.0 and one ion.

Theoretical Absorbance: for each adjacent pair of bases (nearest neighbours) a weight is accumulated using a table. For each internal base a weight is subtracted using another table. Separate tables are used for DNA and RNA.

Calculated factor: this is the calculated molecular weight divided by the theoretical absorbance.

Saving & Printing For details of manual saving and printing see the Saving and Printing section.

#### **Measurement Parameters**

Set the required Base Type (DNA/RNA), whether the base is Phosphorylated, the Primer Concentration (pmoles/ml), Buffer Molarity and Counter Ion. Counter Ion has options for Na (sodium), K (potassium), TEA (triethylammonium) and Other, allowing the user to set the required molecular weight (MW) of the counter ion.

Set the Pathlength and Integration Time you require. Base Sequence allows the user to enter the known base sequence triplets using the buttons A, C, G and T/U. To improve readability a comma is added after each triplet.

Bandwidth (Ultrospec 9000 only) and Lamp Mode can be set as you wish. The Sample Seed entered under Sample will the filename of any automatically saved file.

Set the outputs required in your method. For more information see the section Saving and Printing.

## Taking a Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert a cuvette containing the reference solution in the reference cell holder and a cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

Phosphorylated Counter Ion No Na Primer Conc. 1.000 **Tm Calculation** Pathlength Integration Time 10 mm 0.5 Seconds **Base Sequence** TAA, TAC, GAC, TCA,



CTA, TAG, GG

С

A 4 G

т

Del







## 11.2. Protein Applications

The Ultrospec 7000/8000/9000 contains dedicated methods for both colorimetric protein assays and direct UV measurements.

## 11.2.1. BCA, Bradford, Lowry & Biuret Protein Assays

The BCA, Bradford, Lowry and Biuret protein assays are well established spectrophotometric methods for determining the amount of protein in a sample. The exact choice of the assay depends upon the concentration of protein being measured and the detergents/reducing agents used in purification. Detailed protocols are supplied with all assay kits and should be followed closely to ensure accurate results are obtained. An outline of the protein assays offered by the Ultrospec 7000/8000/9000 is provided below:

Bradford method:	Quantifies the binding of the dye Coomassie Brilliant Blue to an unknown protein and compares this binding to that of different, known concentrations of a standard protein at 595 nm. The standard protein is usually bovine serum albumin (BSA).
Biuret method:	Depends on reaction between Cu <sup>2+</sup> ions and amino acid residues in an alkali solution. The resulting copper complex absorbs light at 546 nm.
BCA method:	Depends on reaction between Cu <sup>2+</sup> ions and amino acid residues. In addition, this method combines this reaction with the enhancement of Cu <sup>+</sup> ion detection using bicinchoninic acid (BCA) as a ligand, giving an absorbance maximum at 562 nm. The BCA process is less sensitive to the presence of detergents used to break down cell walls.
Lowry method:	Depends on quantifying the color obtained from the reaction of Folin-Ciocalteu phenol reagent with the Tyrosyl residues of an unknown protein and comparing with those derived from a

## 11.2.2. Determination Of Protein Concentration Using The Bicinchoninic Acid (BCA) Protein Assay

The principle of the bicinchoninic acid (BCA) protein assay relies on the formation of a Cu<sup>2+</sup>-protein complex under alkaline conditions, followed by reduction of the Cu<sup>2+</sup> to Cu<sup>+</sup>. The amount of reduction is proportional to the amount of protein present. BCA forms a purple-blue complex with Cu<sup>+</sup> in alkaline environments, thus providing a basis to monitor the reduction of alkaline Cu<sup>2+</sup> by proteins. The BCA assay can be used to quantify proteins in the concentration range 0.2 to 1.0 mg/ml. It is compatible with many detergents but not compatible with reducing agents such as dithiothreitol above 1 mM.

standard curve of a standard protein at 750 nm (usually BSA).

## **Getting Started**

It is always advisable to prepare the standard in the same buffer as the sample to minimize any interference effects. BCA assays are routinely performed at 37°C. Color development begins immediately and can be accelerated by incubation at higher temperatures. Higher temperatures and/or longer incubation times can be used for increased sensitivity.

## **Materials Required**

Bicinchoninic Acid Kit for Protein Determination.

Suitable tubes with caps to hold and mix 2.1 ml samples and to heat at up to 60°C.

Plastic disposable cuvettes.

Standard protein solution of known concentration (1 mg/ml).

Incubator or block heater to heat sample tubes.

Preparation Of The BCA Working Reagent.

BCA reagents A and B are available commercially from a number of different sources. Instructions given here are for the kit supplied by Sigma Aldrich, other methods will be similar. Always refer to the manufacturer's instructions.

- 1. Mix 50 parts of Reagent A (a solution containing bicinchoninic acid, sodium carbonate, sodium tartrate and sodium bicarbonate in 0.1 N NaOH, pH 11.25) with 1 part of Reagent B (4% (w/v)  $CuSO_4.5H_2O$ ), preparing sufficient reagent for all the standards and samples. 2 ml of working reagent is required for each sample.
- 2. Mix until the solution is a uniform light green color. The solution is stable for 1 day.

## **Standard Preparation**

- 1. Prepare a series of protein standards ranging in concentration from 0.2 to 1.0 mg/ml such that the final volume for the assay is 0.1 ml. The Ultrospec 7000/8000/9000 can measure up to 9 standards and up to 3 replicates.
- Add 2.0 ml of the BCA working reagent to each standard, vortex gently and incubate using one of the following parameters: 60°C for 15 minutes, 37°C for 30 minutes or room temperature from 2 hours to overnight.
- 3. If required, allow the tubes to cool to room temperature.

## Sample Preparation

- 1. Prepare the unknown samples as described above ensuring that the final volume is 0.1 ml.
- Add 2.0 ml of the BCA working reagent to each sample, vortex gently and incubate using one of the following parameters: 60°C for 15 minutes, 37°C for 30 minutes or room temperature from 2 hours to overnight.
- 3. If required, allow the tubes to cool to room temperature.

## **Creating A Standard Curve**

For BCA measurements *Wavelength* is set to 562.0 nm. *Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* and *Sample* can be set as you wish.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

Set **Calibration** to Standards, *Curve Fit* to Zero Regression and enter *Units* of mg/ml. The number of standards and replicates can be set as you wish (for optimum accuracy it is recommended that the number of standards is  $\geq$ 4 and replicates is >1).

**Note:** With *Calibration* set to Standards the user is required to prepare and measure standards with *Calibration* set to Manual the user inputs both the standard concentrations and absorbances.

Enter the concentration of prepared standards using the numeric keypad.







Set the outputs required in your method. For more information see the section Saving and Printing.

To create the standard curve when using replicates press the replicates button in the bottom right corner to take you to the screen shown below. With *Replicates* off standards can be measured directly as described below.

**Note:** After all replicates have been taken for a standard pressing the replicates icon takes the user to the next standard that was specified in the method.

To create a standard curve using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the first standard/replicate in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the standard measurement.

Continue recording all standards/replicates until the standard curve has been completed. After the standard curve has been collected press the forward arrow to proceed to the sample measurement screen.

To ignore any outlying standard measurements press the tick next in the appropriate row to toggle it to a cross. Any ignored measurement will be automatically removed from the standard curve. These can be reinstated by pressing the cross.

**Note:** Pressing the save method icon before any standards have been measured will save the method parameters only. Recalling a method containing method parameters only will require the user to construct a standard curve before measuring samples.











#### **Taking A Measurement**

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

To view the standard curve whilst on the sample measurement screen, simply press the 'View Curve' icon that appears under the options menu.

**Note:** Saving the method using the Save Method icon that appears on the options menu will save both the method parameters and the standard curve. Recalling a method containing method parameters and standard curve allows the user to measure samples directly.

#### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.

## 11.2.3. Determination Of Protein Concentration Using Direct UV Methods

The direct UV method of protein determination has a number of advantages over traditional colorimetric assays in that it does not rely on an external protein standard and the sample is not consumed in the assay. However the presence of nucleic acid in the protein solution can have a significant effect due to strong nucleotide absorbance at 280 nm. This can be compensated by measuring A260 and applying the equation of Warburg and Christian for the protein crystalline yeast enolase (Equation 1).

Protein concentration (mg/ml) =  $(1.55 \times Abs280) - (0.76 \times Abs260)$  1

Protein concentration = (Factor 1 × Abs280) - (Factor 2 × Abs260) 2

Ultrospec 7000/8000/9000 use default A260 and A280 factors of 0.76 and 1.55, respectively. These factors can be edited so that the equation can be applied to other proteins (Equation 2). Compensation for background, dilution and pathlength can also be entered.

To customize Equation 2 for a particular protein, the A260 and A280 values should be determined at known protein concentrations to generate simple simultaneous equations, which, when solved provides the two coefficients. In cases where Factor 2 is found to be negative, it should be set to zero since it means there is no contribution to the protein concentration due to absorbance at 260 nm.

The A260/A280 ratio also gives an indication of protein purity; a ratio of 0.57 can be expected for pure protein samples.

#### **Background Correction**

- To compensate for the effects of background absorbance caused by turbidity, high absorbance buffer solutions and the use of reduced aperture cells the Ultrospec 7000/8000/9000 can use background correction at 320 nm.
- When used A320 is subtracted from A260 and A280 prior to use so that:

Protein concentration = [Factor 1 ×(Abs 280 - Abs 320)] - [Factor 2 ×(Abs 260 - Abs 320)] Ratio = (Abs 260 - Abs 320) / (Abs 280 - Abs 320)

• The use of background correction can remove variability due to handling effects of low volume disposable cells.



## 11.2.4. Protein UV

## **Measurement Parameters**

Set *Pathlength* and *Dilution Factor* to the required values. Check above to see if background correction is required. The default values for *A260 Factor* and *A280 Factor* are 0.76 and 1.55, respectively; these can be edited by pressing the appropriate box. Set *Units* to encompass the expected concentration of your samples.

You may advance to the next screen at any time by pressing the forward arrow and return to the previous screen by pressing the back arrow.

*Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you wish. The *Sample Seed* entered under *Sample* will the filename of any saved file.

Set the outputs required in your method. For more information see the section Saving and Printing.

## Taking A Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

If *Background* is set to On the A320 result will be included in the left hand column and automatically subtracted from the displayed A260, A280 and A260/A280 results.

#### Saving & Printing

For details of manual saving and printing see the Saving and Printing section.









## 11.2.5. Protein A280

## **Measurement Parameters**

Select the *Mode* you require, Ultrospec 7000/8000/9000 have options for Christian Warburg, BSA, IgG, Lysozyme, Molar Extinction, Mass Extinction and E1% (for Molar Extinction the user will also be required to enter the molar extinction coefficient, and molecular weight and atomic units for Mass Extinction).

Set *Pathlength* and *Dilution* Factor to the required values. Check above to see if background correction is required.

Set *Units* to encompass the expected concentration of your samples. *Bandwidth* (Ultrospec 9000 only), *Integration Time* and *Lamp Mode* can be set as you wish. The *Sample Seed* entered under *Sample* will the filename of any saved file.

Set the outputs required in your method. For more information see the section Saving and Printing.

## Taking A Measurement

To take a measurement using dual beam instruments (Ultrospec 7000/8000/9000) insert the cuvette containing the reference solution in the reference cell holder and the cuvette containing the sample in the sample cell holder and press the take measurement button. The reference measurement will be automatically subtracted from the sample measurement.

If *Background* is set to On the A320 result will be included in the left hand column and automatically subtracted from the displayed A260, A280 and A260/A280 results.

## Saving & Printing

For details of manual saving and printing see the Saving and Printing section.









# 12. SAVING & PRINTING

Ultrospec 7000/8000/9000 UV-Visible spectrophotometers allow users to save and print sample data. This can either be included automatically as a method parameter or performed manually from the sample measurement screen. This section details the saving and printing options offered by all Ultrospec 7000/8000/9000 spectrophotometers.

## 12.1. Saving Sample Data

The Ultrospec 7000/8000/9000 allows users to save sample data in three different formats:

## 12.1.1. Internal

The sample data is saved to the instrument's internal memory format. See the Sample Manager section for details on saving and recalling data from the internal memory.

**Note:** To ensure the optimum performance of the Ultrospec 7000/8000/9000 it is recommended that unwanted data be deleted from the instrument's internal memory at regular intervals.

## 12.1.2. USB

The sample data is saved to a USB memory stick in format that can be read by Ultrospec 7000/8000/9000 spectrophotometers only. Files in this format cannot be opened by Microsoft Excel™ or other spreadsheet programs.

**Note:** Sample data will be saved to the Ultrospec 7000/8000/9000 Samples directory on the USB memory stick; if this directory is not present it will be created.

## 12.1.3. USB CSV

The data is saved to a USB memory stick in comma separated variable (CSV) format allowing it to be opened directly using Microsoft Excel or other spreadsheet programs. Files in this format cannot be opened using the instrument.

**Note:** To view the data displayed in the File Created, Date and Time cells in a recognized format this will need formatting as described below:

File Created: Right click in the appropriate cell and select Format Cells from the list, select Custom dd/mm/yyyy h:mm from the list on the right hand side (see below) and select ok.



Date / Time:

e: Right click in the appropriate cell and select Format Cells from the list, under Category select Date or Time and the desired format from the list on the right hand side (see below) and select OK.



## 12.2. Automatic Saving

The option to save sample data automatically is set under method parameters. With Auto Save set to on, the save location can be set to Internal, USB or USB CSV (the USB options are only available if a USB memory stick is inserted).

The filename given to an automatically saved file will be either the Sample Seed entered in the method parameters or, if the user chooses not to enter a Sample Seed, Default. The Ultrospec 7000/8000/9000 will only save one 'Default' file per application; subsequent saves of files without a sample seed will overwrite the previous default file.

## As sample data is saved when the user exits the application, removing a USB memory stick before exiting will result in loss of data.

**Note:** All files are appended with a unique time and date stamp it is therefore possible to create two or more files sharing the same name.

**Note:** With Overlays  $\geq 2$  in Wavescan and Number of Samples  $\geq 2$  in Kinetics the overlaid data will always be saved automatically to the instrument's internal memory.

## 12.3. Manual Saving

If a method does not require sample data to be saved each time a measurement is taken it is possible to manually save sample data in one of the formats outlined above. This procedure is described below:

After collecting all required sample measurements select save sample data from the options menu on the sample measurement screen to display the dialogue box shown left. The save location and filename are set by pressing the Save to and Sample Name boxes, respectively. If no Sample Name is entered the file will be titled Default.

Note: Any sample data saved manually will override the auto save function

## 12.4. Exporting Data

The Ultrospec 7000/8000/9000 allow users to recall saved sample data from the internal memory or a USB memory stick and save this in another format. This is done as follows:

Recall saved data using Sample Manager and press the save sample button on the options menu to display the save sample dialogue box. Set the desired save location and filename using the Save to and Sample Name boxes, respectively and press the tick to confirm the data export.





# 13. SAMPLE MANAGER

Sample Manager is the application used by the Ultrospec 7000/8000/9000 for saving and recalling data from both the instrument's internal memory and the instrument's USB format. Sample Manager can be accessed from either the main screen using the Sample Manager button (below left) or from within an application using the Sample Manager icon on the options menu (below right).





To recall a saved file, highlight the desired file and press the Sample Manager icon in the right hand corner of the screen. Data will be displayed on the Sample Measurement screen (see Recalled Files below).

With a USB memory stick inserted it is possible to toggle between the internal and USB memories using the icon in the left hand corner of the screen. The location of the displayed data will be indicated by the icon in the top right hand corner.

Sample Manager has been designed to make finding saved files as simple as possible. Therefore it is possible to arrange files alphabetically, by application or by date/time saved by pressing the column headers Sample ID, Application and Date, respectively.

If there are too many saved data files to fit on a single screen this is signified at the top of the screen, e.g. Page 1 of 2. Scrolling through the screens is achieved using the up and down arrows at the bottom of the screen.

**Note:** Sample Manager can only display the first 100 sample data files; if the internal memory or USB stick contains >100 files these can be viewed be deleting or moving (USB only) unwanted data.

## 13.1. Deleting Data From The Internal Memory

To ensure that the internal memory of the instrument does not contain too many unwanted data files, Sample Manager allows you to delete files. This can be done in one of three ways:

Deleting a single file: Deleting multiple files: Highlight the file for deletion and press the delete button. Highlight multiple files and press the delete button.

Deleting all files:

Press the 'Delete all' icon at the bottom of the screen.

**Note:** It is only possible to delete data using Sample Manager accessed from the main screen. Sample Manager does not allow a user to delete sample data files from a USB memory stick; this must be done using a PC.

Sample Manager allows the user to lock files saved to the internal memory to prevent the accidental deletion of files containing precious data. To lock files, highlight the required data and press the lock icon at the bottom of the screen. Locked files are signified by the lock icon in the right hand column. Once a file is locked selecting 'Delete' or 'Delete All' does not clear this data from the instruments' internal memory.

Sample Manager - page 1 of 2 🛛 🔦				
Sample Id	Application	Date	•	
A280	Protein A280	2010/01/26 13:31:05		
Abs 450	Single Wave	2010/01/26 13:21:53		
BCA	BCA	2010/01/26 13:28:23		
Bradford	Bradford	2010/01/26 13:29:11		
Cydye	CyDye DNA	2010/01/26 13:27:05		
DNA	DNA	2010/01/26 13:25:06		
Holmium4	Wavescan	2010/01/26 13:23:19		
Lowry	Lowry	2010/01/26 13:29:54		
<u>_</u> â*	<b>î ≜</b>	Х ≙ ◆		

To unlock a file, highlight the appropriate locked data and press the lock icon at the bottom of the screen.

Note: As sample data files saved to a USB memory stick must be deleted using a PC it is not possible to lock USB sample data using Sample Manager.

To exit from Sample Manager press the 'Exit' icon.

## 13.2. Accessing Sample Manager From The Main Screen

Date

2010/01/26 13:21:53

2010/01/26 13:28:23

2010/01/26 13:29:11

2010/01/26 13:25:06

2010/01/26 13:23:19

2010/01/26 13:29:54

🔒 🔶 🖾

2010/01/26 13:27:05

BCA

Bradford

CvDve DNA

DNA

Wavescan

Lowry

🕯 🔶 🗙 🛛

Ω.

Δ.

Δ

•



When accessed from the main screen Sample Manager will display all files held on the internal memory or USB memory stick and allows the user to lock and delete files saved to the internal memory. As Sample Manager accessed from the main screen allows users to delete files from the instrument's internal memory this option is disabled for Limited users.

To allow post scan manipulation of saved wavescan and kinetics data, files loaded from Sample Manager are loaded directly into Trace Manager (see Trace Manager section for details).

Note: With large numbers of files held on the internal memory there may be a short delay before Sample Manager opens.

## 13.3. Accessing Sample Manager From Within An Application



When accessed from within an application Sample Manager will only display files belonging to that specific application and only allows users to recall saved data and lock files.

## 13.4. Recalled Files

When recalled, sample data files will display the first sample recorded in a measurement. To access all sample data within a recalled file press the Sample box to display a list of all samples contained within the file. Pressing on the desired file will populate the boxes on the sample measurement screen with the saved data. Choosing to measure another sample with an old sample's data displayed simply updates the sample measurement screen and the appropriate boxes.

**Note:** Wavescan and kinetics sample data is recalled using Trace Manager. For details see the Trace Manager section for details.



## below.

described below:

Pressing the Folder box produces a list of available save locations. USB will only appear on the list if a USB memory stick is inserted.

Pressing the Method Name box allows the user to set the desired method name using alpha numeric text entry.

Press the tick to save and exit or the cross to exit without saving.

14. SAVING METHOD

The Ultrospec 7000/8000/9000 allow users to store methods to both instruments internal memory and to USB memory sticks. The procedure for saving methods is

## 14.1. Methods Saved To The Internal Memory

Methods saved to the instrument's internal memory are stored in either the Methods or Favorites folders, both of which are accessed via the main screen. The Biochrom Ultrospec are capable of storing up to 90 methods on the instrument's internal memory.

## 14.2. Methods Folder

The Methods folder is made up of 9 folders each capable of storing up to 9 methods. The method folder icons have been designed to give the user an indication of the number of methods that are stored in that folder.









## 14.3. Renaming Method Folders

14.4. Locking Saved Methods

Using the Rename Folder icon on the options menu (the left hand icon) it is possible to rename any of the 9 method folders using alphanumeric text entry.

Using the lock folder icon on the options menu (the center icon) it is possible to

add pass code protection to any of the method folders. Locked folders cannot be

Within a methods folder it is possible to add pass codes to individual methods and

to lock them from deletion using the lock icon on the options menu (the center

icon). Locked methods can be unlocked by using the unlock icon on the options

menu (the right hand icon), selecting the desired method and entering the correct

#### Methods [ L 8 Methods 1 Methods 2 Methods 3 Δ Δ Δ Methods 4 Methods 5 Methods 6 Options L Δ Methods 7 Methods 9 ♠





Methods - Methods 1

λ

Single Wave

1 Standard Curve

> Options  $\cap$

> > ♠

Ô

2

M

Wavescan

λ

Single Wavelen

L

Kinetics





14.5. Deleting Saved Methods. Within a methods folder it is possible to delete saved methods using the delete icon on the options menu (the left hand icon) and selecting the desired file.

Note: Locked methods must be unlocked to allow deletion.

pass code.

renamed and are indicated by a padlock.

## 14.6. Backing Up Method Folders To USB

With a USB stick inserted it is possible to use the USB icon on the options menu (the right hand icon) to:

Backup Folder:	Copies all methods from a specified folder to a USB memory stick.
Restore Folder:	Copies a backed up method folder from the USB memory stick to the internal memory.
Backup All Folders:	Copies all method folders from the internal memory to a USB memory stick.
Restore All Folders:	Copies all backed up method folders from the USB memory stick to the internal memory.

## 14.7. Favourites Folder

The Favourites folder is capable of storing up to 9 user defined methods. Methods stored in the Favourites folder can be locked and deleted as described above.

## 14.8. Saving Methods To USB

To save files to a USB memory stick follow the procedure described above and select USB in the Folder box. The USB option will only appear in the list if a USB memory stick is inserted.

Methods saved to a USB memory stick will appear in the Ultrospec 7000/8000/9000 Methods folder in the root directory.

**Note:** Although it is possible to save an unlimited number of method files to a USB memory stick only 9 can be displayed on the instrument at any one time. As these will only be read from the Ultrospec 7000/8000/9000 Methods folder, additional files can be stored in other locations









# 15. PRINTING

## Printing Sample Data

The Ultrospec 7000/8000/9000 allows users to print sample data in one of three ways:

Note: Only available printers will be shown in the Print to ... options box.

## 15.1. Internal Printer

Fitting of the internal printer is described in the Ultrospec 7000/8000/9000 Accessories Manual at the end of this manual.

## 15.2. Print Via Computer (PVC)

Print via Computer (PVC) is an application running under Windows™ to enable the Ultrospec 7000/8000/9000 to transfer data into a PC environment. From there the data can be printed or saved in a variety of formats, including graphics and text formats or as an Excel file. PVC can store data either to a common directory or be configured to save to independent directories by both file format and connection.

PVC is capable of supporting several instruments simultaneously, limited only by hardware and the speed of the host system and is able to operate via USB and Bluetooth simultaneously.

Installation and operating instructions for PVC can be found on the PVC CDROM.

## 15.3. Bluetooth®

The Bluetooth accessory is a factory fitted option.

## 15.4. Automatic Printing

The option to print sample data automatically is set under method parameters. With Auto Print set to on, the print location can be set to one of the available options.



## 15.5. Manual Printing

If a method does not require sample data to be printed each time a measurement is taken it is possible to manually print sample data. This procedure is described below:

Set the desired print location in Print to... After collecting all required sample measurements select the Print icon from the options menu on the sample measurement screen.



# 16. Cell Changer

## 16.1. Operation

With a cell changer connected, the Accessories folder under Settings will contain the Cell Changer icon. This provides the user with the means to set the default configuration for the cell changer accessory.

## Prompt User

- Yes: The Cell Changer parameter screen is displayed as part of the application parameters, allowing the user to set cell changer parameters each time an application is run. The default configuration set here pre-selected.
- No: The Cell Changer parameter screen is not displayed as part of the application parameters. The default settings configured in the Accessories folder are used.

The Prompt feature is designed to save the user time and effort configuring each application for the cell changer and will also allow accessory parameter screens to be skipped on loading applications.

## Use As Single Cell

- Yes: Stops the cell changer accessory moving, meaning that the cell position 1 will remain in the sample beam for each measurement. When enabled, the options related to cell changer are removed from the screen. As this option allows the user to use the cell changer as a standard single cell holder this removes the need to keep swapping accessories.
- No: The cell changer accessory will be used.

## Operation

- Automatic: By pressing the Start Measurement button the unit will collect the number of samples specified automatically.Manual: Bu pressing the Start Measurement button the unit will measure
  - anual: By pressing the Start Measurement button the unit will measure the first sample, display a prompt to start next measurement, measure the second sample, prompt display a prompt to start next measurement etc.

## Number Of Samples

Allows the user to set the number of samples to be measured using the numeric entry box. In Fixed Wavelength applications, if the number of samples is >8 the user will be prompted to load the cell changer after the first 8 samples. In Wavescan and Kinetics the maximum number of samples is 8.

**Note:** Number of Samples refers to the number of samples only and does not include standard or reference measurements.

## Prompt For Cell Pos

- On: After pressing the Start Measurement button the software will display the positions each sample should be placed.
- Off: The unit will perform the measurement without displaying this message.

The Cell Changer parameter screen that is displayed in the application method parameter screens is shown right (with use as single cell on or off). All options are as described above.

Cell Changer		
Prompt User	Number of samples	
Yes	1	
Use as single cell	Ref after sample	
No	No	
Operation	Prompt for cell pos	
Automatic	Off	
×	$\checkmark$	

Use as single cell No Ref after sample No No Operation Automatic Off Number of samples 1 Cell Changer Parameters Use as single cell Yes	Cell C	hanger
No     No       Operation     Prompt for cell pos       Automatic     Off       Number of samples     1       1     Cell Changer Parameters       Use as single cell     Yes	Use as single cell	Ref after sample
Operation     Prompt for cell pos       Automatic     Off       Number of samples     1       1     Cell Changer Parameters       Use as single cell     Yes	No	No
Automatic Off  Number of samples 1  Cell Changer Parameters Use as single cell Yes	Operation	Prompt for cell pos
Number of samples 1 Cell Changer Parameters Use as single cell Yes	Automatic	Off
1 Cell Changer Parameters	Number of samples	
Cell Changer Parameters Use as single cell Yes	1	
Cell Changer Parameters Use as single cell Yes	×	5
Use as single cell Yes	Cell Change	er Parameters
Yes	Use as single cell	
	Yes	
	-	•

### Taking A Measurement With A Cell Changer

When using the cell changer the software will  $\underline{not}$  display the Take Reference button.

Ultrospec 7000/8000/9000 As these instruments continually reference against the reference cell holder there is no requirement to perform a reference measurement before collecting sample data.

The icon in the top right corner of the sample measurement screen indicates the cell position that is in the sample measurement beam.

## 16.2. Saving Methods & Results

Any method saved with the cell changer connected will save the inputted accessory parameters. If a user attempts to use a cell changer method without this accessory connected the software will display the warning, 'The following accessories are missing: Cell Changer The cell changer must be fitted before this method can be loaded.'.

## 16.3. Incorrect Fitting

If the cell changer is fitted incorrectly it will not be recognized during the instruments' start up calibration tests. If this occurs the instrument should be switched off, disconnected from the mains and the accessory refitted, paying close attention to the internal accessory connector.

## 16.4. Using The Cell Changer To Create A Standard Curve

Pressing the take measurement button when on the Replicates screen will re-measure the highlighted replicate only. This measurement will always be made from cell position 1 of the cell changer.

Pressing the take measurement button on the construct standard curve screen will re-measure all standards whether measuring replicates or not.







# 17. Sipper

The Sipper and Flowcell accessories for the Ultrospec range of instruments are designed as an aid to speed up sample handling. The system comprises a peristaltic pump head that is controlled by the software to sip a known, calibrated volume of sample and move this into a Flowcell which is in the sample position of the spectrophotometer.

The Sipper can be set up and used within all applications and Sipper parameter are stored as part of an application method.

## 17.1. Setting Sipper defaults

It is simple to customize the default parameters of the Sipper for a particular installation by setting all of the default values. Every application will then load these defaults each time the Sipper page is opened however they can be changed within the application. These default values are set within the instrument under the Settings, Accessories, Sipper Icons and are stored even when the instrument is switched off.



#### Sip Volume

• Amount of sample to be sipped up. Range of values = 0.3 ml to 50 ml.

**Note:** working with volumes less than 1 ml will require careful setting of the air gap to minimize cross contamination between samples.

#### Air-gap

 Gap between successive samples which is set by pumping air after the tube has been removed from the sample liquid – set in cm of tubing. Range of values = 0 to 200 cm.

The Sipper has the ability to program an air gap between samples to help minimize cross contamination and to allow users flexibility to work with different volumes and lengths of Sipper inlet tubing. When the start button is pressed the Sipper pumps up the volume of sample that is set within the method and the user is then prompted to remove the tubing from the sample. After a short delay the pump starts again and the sample is moved along the tubing and air is sucked into the end of the tubing. The default air gap set in the software is 40 cm which has been set to work with a 65 cm length of tubing on the Flowcell inlet. If a different length of tubing is used then a different default air gap will need to be set. To calculate this default air gap:

- Measure the length of tubing connected to the Flowcell tubing inlet and take 20 cm from this value.
- From the home page go into a single wavelength method and enter through until you get to the Sipper parameter page. Set the volume to 1 ml and the air gap to the value measured above. Set the settle time to 1 second and Auto read to off.
- Sip a sample of water, remove the tube from the sample when prompted and note the position of the end of the sample slug – you need this to be stopping above the inlet position to the Flowcell. For a 1 ml sample volume you should have this stopping 15-25 cm above the Flowcell inlet. To move the position of the sample press the menu icon, parameters icon, step through the method until you reach the Sipper parameters and adjust the air gap value. Increasing the value will move the end of the sample closer to the Flowcell inlet, lowering the value will move it further away.
- Once the value has been determined exit the Single wavelength application, go to Settings, accessories, Sipper and enter this value in the default air gap box.

#### Settle time

• Amount of time before the system is ready to measure after the sample has entered the Flowcell. Value entered in seconds. Range of values = 1 to 30 seconds

### Sip Mode

 Choice of Waste or Return. Waste mode turns the pump head in one direction and the sample goes to waste after measurement. If return is set then pressing the start button a second time will reverse the motor and the sample will be pumped back out of the inlet tube into the sample vessel. Cross contamination between samples will be worse if return mode is used.

#### Auto Read

• Auto read on will automatically start the measurement in the current application after the settle time has passed. If Auto read is off the user will manually have to press the reference or run icon for each measurement.

If you press the Settings, accessories, Sipper icons as well as the defaults that have been set above there are other icons at the bottom of the screen for Calibration of the Sipper volume and washing of the Sipper.

## 17.2. Calibration

The Sipper software contains a routine that allows users to calibrate the Sipper to ensure that the volume of liquid set within the method matches the actual volume being sipped. Before starting the calibration routine check that it is actually necessary:

- Sample volume set up a single wavelength method which includes the Sipper and set the sip volume to 1 ml. Fill a 10 ml measuring cylinder, sip 5 times from the measuring cylinder taking notice of the messages regarding removing the tube for the air gap. After 5 sips divide the volume of sample sipped by 5 – the value should be approximately 1 ml. If the value varies by more than 0.2 mls then re-calibrate the Sipper as below:
- Fill a 10 ml graduated measuring cylinder with water up to the 10 ml line. Press the calibrate icon, put the end of the sample tube into the measuring cylinder, press the start button and the Sipper will start pumping. Once this has finished repeat 4 more times. Read off from the measuring cylinder the actual amount that has been sipped and enter it into the box on the screen and then press the Tick icon to exit.

If you need to rest the calibrations to the default value then sip once and enter the value of 1 ml.

## 17.3. Wash mode

This enables you to wash the Sipper at the end of a batch of samples. Press the wash icon, dip the sample tube into water and allow it to flush the system.

#### Using the Sipper within an application

All Sipper parameters can be set as part of a method, the following screen will appear as part of the application method when the Sipper is connected. The screen will open using the default Sipper values stored under the Settings, Accessories, Sipper Icons.

The process of using the Sipper is as follows:

- 1. Choose the application that you want to run by pressing the relevant icon.
- 2. Set up the instrument parameters, data handling parameters and then Sipper parameters. The Sipper parameters shown will be the ones that are set on the Sipper default page. These Sipper parameters can be changed as part of this application method and the new values saved as part of the method.
- **3.** Put the sample tubing into the reference solution and press the silver button on the front of the Sipper. System will sip the set volume and then prompt you to remove the sample tubing.
- **4.** System will then pump the set air gap and then wait the set settle time. If auto read is on the system will automatically reference, if it is off then press the clear cuvette icon to reference.





- 5. Put the sample tube into the sample, press the silver button on the front of the Sipper. System will sip the set volume and then prompt you to remove the sample tubing.
- **6.** System will then pump the set air gap and then wait the set settle time. If auto read is on, the system will automatically run the sample, if it is off then press the orange cuvette icon to run the sample. Repeat for subsequent samples.

## 17.4. Cleaning/Maintenance/Troubleshooting

To ensure good performance from your Sipper system it is essential that routine maintenance is carried out.

### After batch of samples

• Run the Sipper Wash mode pumping deionized water, for 1 minute then remove sample tubing and pump air until all liquid is pumped into the waste vessel.

#### At end of day

• Open up pump head to relax pressure on pump head tubing.

## Weekly

- Check pump head tubing for signs of wear replace if necessary.
- Check all connections for leaks.
- Depending upon types of sample being run it may be advantageous to clean the Flowcell properly (e.g. if "sticky" protein solutions are being run). Unscrew the 2 connectors, take the Flowcell out of the instrument, fill with a detergent solution and leave to soak overnight. Empty and then flush with deionized water for 1 minute using the wash mode.

Problems with Sipper systems are generally caused by wrongly set parameters or poor maintenance. Typical faults and their cures include:

### Sample not being sipped

- Check pump head closed.
- Check pump head tubing for damage.
- Check Flowcell is connected and connectors are tightened.
- Check PTFE tubing is pushed firmly into pump head tubing.
- Check PTFE tubing is not crushed or damaged.
- Check Flowcell for blockages.
- Check waste pipe for blockages.

## Sample being pumped right through Flowcell

• Check air gap and reset to correct value.

#### **Bubbles in Flowcell**

- Check Flowcell connectors are tightened.
- Clean Flowcell with detergent solution.

#### Inconsistent readings

- Check volume being sipped try increasing the volume.
- Adjust air gap so the end of the sample is closer to the Flowcell inlet more of the sample flushes the previous sample out of the Flowcell.

### Wrong volume being sipped

• Check Sipper calibration.

# 18. TECHNICAL SPECIFICATIONS

Parameter	Ultrospec 7000 Specification	Ultrospec 8000 Specification	Ultrospec 9000 specification
Configuration	Dual Beam	Dual Beam	Dual Beam
Lamp	Xenon	Tungsten/Deuterium	Tungsten/Deuterium
Wavelength Range	190 nm to 1100 nm	190 nm to 1100 nm	190 nm to 1100 nm
Wavelength Accuracy	+/- 0.5 nm across wavelength range	+/- 0.3 nm across wavelength range	+/- 0.3 nm across wavelength range
Wavelength Reproducibility	+/- 0.1 nm	+/- 0.1 nm	+/- 0.1 nm
Bandwidth	2 nm	1 nm	Variable (0.5 nm, 1 nm, 2 nm, 4 nm)
Toluene in Hexane EP Resolution		>2.0	>2.0 at 1 nm bandwidth
Stray Light	<0.050% T at 220 nm using Nal and 340 nm using NaNO $_3$	<0.025% T at 220 nm using Nal and 340 nm using NaNO <sub>3</sub> <1% T at 198 nm using KCl	<0.025% T at 220 nm using Nal and 340 nm using NaNO <sub>3</sub> <1% T at 198 nm using KCl
Photometric Range	-4 A to 4 A	-4 A to 4 A	-4 A to 4 A
Photometric Accuracy	+/-0.002 A at 0.5 A +/-0.004 A at 1 A +/-0.006 A at 2 A	+/-0.002 A at 0.5 A +/-0.004 A at 1 A +/-0.006 A at 2 A	+/-0.002 A at 0.5 A +/-0.004 A at 1 A +/-0.006 A at 2 A
Photometric Reproducibility	+/-0.002 A at 1A	+/-0.002 A at 1A	+/-0.002 A at 1A
Scan Speed	>2400 nm/min	>2400 nm/min	>2400 nm/min
Zero Stability	+/-0.001 A/hr	+/-0.0003 A/hr in precision mode	+/-0.0003 A/hr in precision mode
Noise	< 0.00020 A RMS @ 0A at 700 nm and 500 nm over 20 measurements	< 0.00005 A RMS @ 0A at 700 nm and 500 nm over 20 measurements	< 0.00005 A RMS @ 0A at 700 nm and 500 nm over 20 measurements

# 19. CHANGING TUNGSTEN AND DEUTERIUM LAMPS



Lamps should only be removed after they have cooled completely, attempting to remove lamps before they have cooled may result in injury.

**1.** Switch off the Ultrospec 8000/9000 and ensure that the instrument is disconnected from the mains.

**2.** From the back of instrument remove the screws securing the lamp chamber (this is located above the fan grating).

**3.** Remove the lamp chamber lid to reveal both the Tungsten and Deuterium lamps (see below).







Tungsten Lamp

Deuterium Lamp



LAMP CHAMBER OF THE Ultrospec 8000 & 9000

## Removing & Replacing The Tungsten Lamp

To remove the tungsten lamp, pull the metal clip away from the tip of the lamp and carefully pull the lamp to remove.

To insert the new lamp, pull the metal clip away from the housing to provide space for the new lamp to be inserted and carefully insert the pins of the lamp into the lamp housing. When inserted in the housing carefully place the metal clip on the tip of the lamp.

**Note:** When replacing a tungsten lamp, the lamp envelope must not be touched. If the envelope is touched during installation the lamp must be thoroughly cleaned with a light cloth.

### Removing & Replacing The Deuterium Lamp

The Deuterium lamp is removed by carefully pulling the lamp upwards. To insert the new lamp, line up the lamp pins with the holes in the lamp housing and push into place. The Deuterium lamp can only be inserted in the correct position.

**Note:** After installation the Deuterium lamp must be thoroughly cleaned with a light cloth to remove any marks.

#### Replacing The Lamp Chamber Lid

After the new lamp has been installed the lamp chamber lid should be replaced and screwed in place. The instrument will now be ready to use.

**Note:** A new baseline should always be measured after installing a lamp. This procedure is described in "8.6. Instrument Settings" on page 18. The baseline should only be collected after the lamp has warmed up sufficiently (at least 30 minutes).

## **Resetting Lamp Life**

With new lamps installed it is necessary to reset the lamp life on the software, this is done as follows:

From the main page, select Settings followed by Instrument Settings.



Lamp Settings			
	Deuterium	Tungsten	
Lamp State	On	On	
Installed Date Installed Time	1//1//1980 00:00:00	1//1//1980 00:00:00	
Original Energy Current Energy	0.0 0.4	0.0 0.0	
Turned On Hours On Total Hours On	1 0 0	2 0 0	
	<b></b> <sup>0</sup>	Q®	
1 i ·	2		×

Select the lamp status icon at the bottom of the screen and press the appropriate Reset Lamp Life icon to reset the lamp life information for the desired lamp. The icons are detailed in the Table of Icons section.

**Note:** Only users with Administrator privileges are capable of resetting lamp life information. For PC variants, lamp life can be reset using Datrys software. See Datrys user manual or on-line help for more details.

# 20. TABLE OF ICONS

Icon	Title	Function
♦	Back arrow	Returns the user to the previous screen.
•	Forward arrow	Advances the next screen in a sequence.
t	Open options menu	Displays the relevant options menu, the exact content of the menu will depend upon the location.
+	Page down	Allows the user to navigate to the next page.
+	Page up	Allows the user to navigate to the previous page.
ŧ	Cursor left	Used in the wavescan and kinetics applications to move the cursor left. The position of the cursor and the corresponding x and y values are displayed above the scan.
1	Cursor right	Used in the wavescan and kinetics applications to move cursor right. The position of the cursor and the corresponding x and y values are displayed above the scan.
0	Toggle data viewed/reset	When used in the CyDye DNA application this toggles the data displayed between DNA and dye data. Under Instrument Settings this will reset a value.
* Ē	Delete all	Deletes all saved data/methods. Delete all is a two stage process.
Â	Delete	Deletes the highlighted data/method. Delete is a two stage process.
	Thresholds	Used in Equation Editor. This allows users to add thresholds (pass/fail) limits to their results.
$\checkmark$	OK/accept	Used to confirm/accept any changes.
Х	Exit	Exits from an application or screen. If exit is selected without saving, any changes will be lost.
	Lock	Locks sample data file/method folder against accidental deletion.
$\sim$	Unlock	Unlocks a locked method folder.
_	Instrument	Used in Sample Manager to access sample data files stored on the instrument's internal memory or to indicate that the data being displayed is from the internal memory.
6	Add linear section	Used in the kinetics applications to view a line of best fit on the scan.

lcon	Title	Function
to	Section	Used in kinetics applications to view data in a specific section. It is possible to add t0 through to t7.
	Method folder	Methods saved internally can be stored in one of nine method folders (each capable of storing up to nine individual methods).
-	Rename method folder	Accessed via the options menu on the methods screen, this allows the method folder to be renamed.
٨A	Toggle view scan on or off	Used in the nucleic acid applications DNA, RNA and Oligo this allows the user to view a survey scan of the last sample run (in the region 220–320 nm).
	Toggle auto print on or off	Accessed via the options button on the sample screen. Pressing this button toggles auto print on and off.
	View method parameters	Accessed via the options button on the sample measurement screen. Pressing this button takes the user back to the method parameters.
۲. ۲.	Print data	Accessed via the options button on the sample measurement screen. Pressing this button prints the sample data.
	Save method parameters	Accessed via the options button on the sample measurement screen. This allows a method to be saved to a location specified by the user.
	Save sample data	Accessed via the options button on the sample measurement screen. This allows sample data to be saved to a location specified by the user.
	Sample Manager	Accessed via the options button on the sample measurement screen. Displays the sample data held in either the internal memory or on a USB memory stick.
i	Instrument Information	Accessed via instrument settings. This displays instrument information such as product name, serial number etc.
۶	Instrument Settings	Accessed via instrument settings. This allows the user to set the default bandwidth, save new baselines and view the date of the last service.
	Instrument Status	Accessed via instrument settings. This displays the current status of the instrument.
0	Lamp Settings	Accessed via instrument settings. This displays the lamp status and allows the user reset the lamp life (8000 & 9000 only).
	Save instrument options	Accessed via instrument settings. This saves the instrument options.
٦	Cell changer number	Displayed on the status bar. This indicates that the cell changer is connected and the cell position that is in the sample beam.
$\ast$	Auto print - Bluetooth	Displayed on the status bar. This indicates that the instrument will automatically print all sample data to the Bluetooth device.
Ð	Auto print - internal printer	Displayed on the status bar. This indicates that the instrument will automatically print all sample data to the internal printer.

lcon	Title	Function
÷	Auto print - USB	Displayed on the status bar. This indicates that the instrument will automatically print all sample data to the USB printer.
Ô	Deuterium lamp off	Displayed on the status bar. This indicates that the Deuterium lamp is off.
M	Deuterium lamp fail	Displayed on the status bar. This indicates that the Deuterium lamp has failed.
Ê	Deuterium lamp on	Displayed on the status bar. This indicates that the Deuterium lamp is on.
Î	Deuterium lamp warming	Displayed on the status bar. This indicates that the Deuterium lamp is warming up.
X	Tungsten lamp fail.	Displayed on the status bar. This indicates that the Tungsten lamp has failed.
Q	Tungsten lamp off	Displayed on the status bar. This indicates that the Tungsten lamp is off.
Q	Tungsten lamp on	Displayed on the status bar. This indicates that the Tungsten lamp is on.
Â	Tungsten lamp warming	Displayed on the status bar. This indicates that the Tungsten lamp is warming up.
$\mathbf{\widetilde{M}}$	Xenon lamp failed	Displayed on the status bar. This indicates that the Xenon lamp has failed.
⊠	Instrument busy	Displayed on the status bar. This indicates that the instrument is performing a measurement.
*	Printing to Bluetooth device	Displayed on the status bar. This indicates that the instrument is printing to a Bluetooth device.
B	Printing to internal printer	Displayed on the status bar. This indicates that the instrument is printing to the internal printer.
÷	Printing to USB	Displayed on the status bar. This indicates that the instrument is printing to a USB printer.
1	Toggle view standard curve	When used in standard curve applications i.e. Lowry protein assay, this toggles the sample measurement screen to display or hide the standard curve.
(0)	Replicates	Used in standard curve applications to collect data from a group of replicates.
7	Take sample measurement	Commences a sample measurement.
Ø	Take reference measurement	Commences a reference measurement.

lcon	Title	Function
◄	Backspace/delete	Used in text entry mode to move the cursor backwards (left) and delete any unwanted characters.
►	Forward space	Used in text entry mode to move the cursor forwards (right).
abo	Lower case	Selects lower case letters when in text entry mode.
λμ!	Symbols	Selects symbols when in text entry mode.
ABC	Upper case	Selects upper case letters when in text entry mode.
123	Numeric	Selects numeric entry when in text entry mode.
1	Trace Manager	Used in the wavescan and kinetics applications. This allows the user to overlay up to 8 samples data files and to choose what type of data is displayed i.e. raw data, smoothed data, 1st derivative.
<b>~</b>	USB	Used on the Sample Manager screen to access data stored on a USB memory stick. Used in the options menu in the methods folder to allow the user to backup methods to a USB memory stick.
	Add user	Displayed under User Access, this allows anyone with Administrator privileges to add another user to the instrument.
8	Delete user	Displayed under User Access, this allows anyone with Administrator privileges to delete users from the instrument.
8	Edit user	Displayed under User Access, this allows anyone with Administrator privileges to edit currently users' parameters.
Ð	Zoom in	Used in the wavescan and kinetics applications. This allows the user to zoom into a specific region of a scan.
Θ	Zoom out	Used in the wavescan and kinetics applications. This allows the user to zoom out and return to the original scan.
# 21. GLOSSARY OF BOXES

Вох	Function
Auto-Print On	Toggles between on and off. Used in all applications to set whether sample data is printed automatically or not.
Auto Save On	Toggles between on and off. Used in all applications to set whether sample data is saved automatically or not.
Background On	Can be toggled on and off. Used in nucleic acid and protein measurements to subtract the absorbance value at 320 nm. This is done to allow for the effects of turbidity, high absorbance buffer solutions and the use of reduced aperture cells.
Background Wavelength 320.0 nm	Set by numeric entry. Used in CyDye DNA measurements only and enables the user to specify the wavelength of background correction.
Bandwidth 2nm	For variable bandwidth instruments (Ultrospec 9000) this gives the option to set the bandwidth to 4, 2, 1 or 0.5 nm for a measurement. For fixed bandwidths instruments (Ultrospec 7000 & 8000) this box will be greyed out.
Base Sequence       A     C     G     T     Del	Used in T <sub>m</sub> calculation to set the sequence of bases. The bases A, C, G, T can be added in DNA mode and the bases A, C, G, U can be added in RNA mode. Bases are grouped in threes to improve readability.
Base Type DNA	Used in T <sub>m</sub> calculation to toggle between DNA and RNA base types.
Brightness 8	Used in User Interface in Settings to set the brightness of the screen.
Buffer Molarity 0.100	Numeric entry, used in T <sub>m</sub> calculation only. Buffer molarity = buffer molarity + total molarity of salt (moles / L).
Calibration Standards	Used in all standard curve applications to set the method used to collect the standard curve. Options are for Standards or Manual.
Correction Factor 0.060	Set by numeric entry, used in the CyDye DNA application only. This is the correction factor applied to the absorbance of the dye at a specified wavelength.
Counter Ion Other	Used in T <sub>m</sub> calculation only this allows the user to add the type of counter ion used. The options are sodium (Na), potassium (K), triethylammonium (TEA) or other. If other is selected the molecular weight of this counter ion must be added using Other MW.
Curve Fit Regression	Used in all standard curve applications this is the curve fit that will be applied to the standards' absorbance values. Options are Zero Regression, Regression, Interpolation and Cubic Spline.
Custom Dye Name	Alphanumeric entry for custom dye name, used in CyDye DNA only.
Delay 0	Numeric entry. This is the delay required before a kinetic measurement commences.
Dilution Factor	Numeric entry. This is used in nucleic acid and protein applications to compensate for the absorbance of highly concentrated samples.

Вох	Function			
Draw Peaks On	Used in wavescan only, this switches display of peak cursors on and off. Peak cursors show vertical dashed lines displaying the measured peak height and horizontal dashed lines showing the peak width.			
Duration 1:30	Numeric entry. This is the duration over which a kinetic measurement is performed.			
Dye 1 Name Cy3	Used in CyDye DNA application only. This is the first dye type used in the measurement.			
Dye 2 Name Alexa Fluor 546	Used in CyDye DNA application only. This is the second dye type used in the measurement (where applicable).			
Extinction Coefficient 150.0 E+3	Numeric entry for CyDye DNA application only. This is the extinction coefficient of the specified dye. Note for dyes included in the Ultrospec software these values are not editable and the box will be greyed out.			
Factor 1000.0	Numeric entry to 3 decimal places. In concentration and nucleic acid measurements multiplying absorbance readings by this factor gives the concentration value. In kinetic measurements the result is calculated by multiplying this factor by absorbance, delta absorbance or the slope.			
Feature Detection Sensitive	Used in wavescan measurements only. This determines the sensitivity of the peak or valley detection i.e. sensitive will detect more peaks or valleys than course. Options are; Off, Coarse, Sensitive or Custom (when custom is selected the minimum peak height and width must be entered).			
Feature Sort Wavelength	Foggles between wavelength and absorbance, used in wavescan measurements only. This determines how features will be ordered in the peak/valley table below the scan.			
Feature Type Peaks	Toggles between peaks and valleys, used in wavescan measurements only. This determines what feature type will be detected.			
Group Limited	Used in User Access to set the group a user will belong to and what access they will be granted.			
Integration Time 1 second	Used in all applications. This is the duration the instrument will take a reading at an individual wavelength. The longer the integration time, the greater the signal to noise ratio and the greater the accuracy.			
Interval 1.000	Numeric entry. This is the interval at which serial kinetics readings will be taken.			
Lamp Mode Precision	Used by the Ultrospec 8000 & 9000 in all applications to toggle between Precision and Pulse mode.			
Max Wavelength 500 nm	Numeric entry. This is the upper limit of a wavescan measurement. Note that when using the Ultrospec 7000/8000/9000 the max wavelength must always be greater than the min wavelength by at least the step value.			
Measurement Mode Serial	Toggles between the kinetics measurement modes, serial and parallel.			
Min Wavelength 400 nm	Numeric entry. This is the lower limit of a wavescan measurement. Note that when using the Ultrospec 7000/8000/9000 the max wavelength must always be greater than the min wavelength by at least the step value.			
Mode Absorbance	Used in the Single Wavelength, Kinetics and Protein A280 applications to set the required measurement mode.			
Nucleic Acids dsDNA(260nm)	Used in CyDye DNA application only, this sets the nucleic acid used in the measurement.			

Вох	Function		
Number of Dyes 2	Toggles between 1 and 2. Used in CyDye DNA application only to set the number of dyes used in the measurement.		
0ther MW 1.000	Numeric entry. This option is only used if the counter ion type in the ${\rm T}_{\rm m}$ calculation is set to Other.		
Password 1000	Used in User Access to set a password for new users.		
Pathlength 10 mm	Used in the nucleic acid applications, Protein A280 and Protein UV. This is the pathlength of the cuvette used in the measurement. Options are for 10, 5, 2 and 1 mm.		
Phosphorylated No	Toggles between yes and no. Used in $\rm T_m$ calculation to set if the sample to be measured is phosphorylated or not.		
Primer Conc. 1.000	Numeric entry to 3 decimal places. Used in $\rm T_m$ calculation to sets the primer concentration in pmole/mL.		
Print to Internal Printer	Used in all applications to set the desired print location. Only available printers are displayed.		
Prompt between λ Off	Toggles between On and Off, used in Equation Editor only. With Prompt on the measurement will proceed as follows: measure wavelength 1, prompt for sample, measure wavelength 2, prompt for sample etc.		
Replicates 3	Used in all standard curve applications. This is the number of times a standard measurement is repeated before the mean of these values is plotted on the standard curve. Options are off (1 measurement), 2 or 3.		
Sample Overlays 4	Used in wavescan measurements, this determines how many samples will be overlaid on the graph. Options are off, 2 to 8.		
Samples 1	Used in kinetic measurements, this sets the number of samples that will be measured during the method. Options are 1, 2 or 3.		
Save to USB	Used in all applications to set the desired save location. Only available locations are displayed.		
Scan On	Toggles between on and off. Used in DNA/RNA/Oligo measurements this sets whether, as a further analytical tool, the instrument will also perform a scan over the wavelength range 220-320 nm.		
Scan Speed Medium	Used in wavescan measurements, this determines the speed at which the scan is performed. Options are Slow, Medium, Fast or Integration. Note that when integration is selected the user will be prompted to set the integration time (see box above).		
Screensaver Off	Used in User Interface in Settings to set the time before the screen-saver will be displayed.		
Show Login Yes	Used by the Default Administrator in User Access to set if user login will be displayed or not.		
Standards 6	Used in all standard curve applications. This is the number of standards that will be used to create the standard curve; options are from 1 to 9.		
Std. 1 10.00	Numeric entry to 2 decimal places. Used in all standard curve applications this is the concentration of the standard.		
Step 1.0nm	Used in wavescan measurements, this determines the interval at which wavescan measurements are made (in nm). Options are for 0.1, 0.2, 0.5 or 1 nm. Smaller intervals give greater detail but lead to longer scan times.		

Вох	Function	
Text Entry Mode	Used in User Interface in Settings to set the text entry mode used for alphanumeric text entry.	
Units         Used in applications that give a concentration result. Units are entered via either           alphanumeric entry or from a list of options.		
User Name	Used in User Access when creating a new user.	
Volume (μl) 2.000	Numeric entry to 3 decimal places, this is used in the CyDye DNA application and is the volume of the probe in $\mu\text{L}.$	
Wavelength 450.0 nm	Numeric entry to 1 decimal place and is used in all fixed wavelength applications to determine the wavelength at which the measurement will be performed.	
λ Max 550.0 nm	Numeric entry to 1 decimal place. Used in the CyDye DNA application, this is the wavelength at which the absorbance of the dye will be measured. Note for dyes included in the Ultrospec software these values are not editable and this box will be greyed out.	
<b>Ү Мах</b> 2.000	Numeric entry to 3 decimal places. This is the maximum value of the Y axis shown during a kinetics measurement. Note the graph will automatically re-scale at the end of the measurement to give the optimum Y max.	
Y Min 0.000	Numeric entry to 3 decimal places. This is the minimum value of the Y axis shown during a kinetics measurement. Note the graph will automatically re-scale at the end of the measurement to give the optimum Y min.	

# Ultrospec<sup>™</sup> 7000/8000/9000

# Accessories



# 1. INSTALLATION

Before installing any accessory in the Ultrospec 7000/8000/9000 you must ensure that the instrument is disconnected from the mains.

# 2. PASSIVE ACCESSORIES

Accessories that do not interact with the instrument's software are named passive accessories. Their installation is described below:

To allow easy access to the thumbscrew at the front of the accessory open the cell chamber lid and lift out the front panel (below left).



With the front panel removed the accessory is removed and replaced as outlined below:

1. Loosen the front thumb screw by turning anti-clockwise.

**2.** With the thumb screw fully unscrewed, lift the accessory and gently pull forward.

**3.** To insert the desired accessory, lift the front end, slide into the brackets at the back of the cell chamber (indicated by the two arrows) and tighten the thumb screw by turning clockwise.

With the accessory inserted simply replace front panel, switch on the instrument and it will be ready to use.







# 2.1. Single Cell Holder (supplied with unit)

The single cell holder is designed to take measurements of 10 mm pathlength cuvettes. All Ultrospec 7000/8000/9000 are supplied with a single cell holder as standard.

# 2.2. Variable Pathlength Cell Holder 80-7100-05

The variable pathlength cell holder is designed to use 1, 2, 5, 10, 20, 50 & 100 mm pathlength cells. To accommodate different pathlengths the clips should be moved to secure the cell and spacers used as required.

# 2.3. Water Thermostatted Cell Holder 80-7100-03

The water Thermostatted cell holder is designed to work with water circulators to achieve constant temperatures in the range 5 to 80°C. The accessory has to be connected to an external fluid circulator so that the water flows into the bottom of the cell holder (the right hand inlet) and flows out through the top of the cell holder (the left hand outlet).

# 2.4. Film Holder 80-7100-12

The film holder is designed to measure the absorbance of solid materials including thin flexible sheet materials such as plastic, acrylic, gelatine. Applications also include the quality control of tinted optical lenses.

# 2.5. Test Tube Holder 80-7100-11

The test tube holder allows samples to be measured in a test tube of between 10 and 24 mm diameter and up to 160 mm tall.











# 2.6. Peltier Controlled Fluid Circulation 80-7100-50

# 2.6.1. General Information

## 2.6.1.1. General Warnings

- Check the voltage selector
- Always disconnect the mains plug before starting any work inside the instrument
- Set the temperature to the required values on power on
- If using substances dangerous for health and the environment, observe the laws and the standards in force in the laboratory where this instrument is used
- If using flammable or explosive substances take the necessary precautions

## 2.6.1.2. Quality, Reliability and Safety

This equipment has been designed with emphasis on QUALITY, RELIABILITY and SAFETY, and we will accept responsibility for these aspects only if the following conditions are met:

- Electrical installation of the room or building in which the equipment is to be used must comply with regulations specified in the country where the equipment is to be used.
- The equipment must be used in accordance with the instructions for use provided by Biochrom.
- All modifications and repairs to the equipment must be carried out by Biochrom or its authorized agents.
- Modifications must not be carried out unless they conform to approved Engineering Service Information issued according to the appropriate Biochrom procedure.
- The equipment installation must be carried out in accordance with local requirements for responsibility and warranty.

# 2.6.2. Instrument Description

The Controller is a Peltier effect thermostatting system using an internal sealed liquid circuit allowing full compatibility and connection to all thermostattable accessories.

The Peltier effect is based on the reversed thermoelectric effect. This is a phenomenon in which a temperature difference is caused by a flow of current through a closed circuit consisting of two different metals.

The Controller provides temperatures ranging from 20°C to 60°C with  $\pm$  1°C stability and  $\pm$  0.1°C repeatability. The features of the Controller include:

- Peltier effect for heating and cooling the liquid of circulation
- Internal flow pump to circulate the liquid
- The Controller can operate in either manual or remote control via a USB port
- The Controller is supplied with one set of thermally insulated connection tubes to connect the Controller to any thermostattable accessory
- The Controller is supplied with a water level sensor to ensure correct operation
- The Controller is supplied with high temperature sensor to avoid damaging the system and malfunctions



## 2.6.2.1. Controller front and rear panel

The front and the rear panel have grids to dissipate the heat generated in the unit. Two fans force the air movement. These grids must be free to guarantee the instrument specifications.

The front panel features the display with keypad.

The rear panel features the mains power connection with mains voltage selector, the USB connector for remote control, the mains switch and the "Water Fill" inlet and outlet connections.

#### 2.6.2.2. Specifications

#### 2.6.2.2.1. Keypad/Display Unit

The keypad/display unit has a LCD screen. All programmable parameters can be displayed in specific menus.

Three buttons are used to access the various functions and change the set parameters.

#### 2.6.2.2.2. Programmable Parameters

• Temperature

#### 2.6.2.2.3. Displayed Values

- The current thermostat temperature
- The set temperature
- The heating/cooling power

#### 2.6.2.2.4. Operating Specifications

- Temperature range from 20°C to 60°C
- Temperature accuracy better than  $\pm$  0.15°C from 20°C to 60°C
- Temperature repeatability typically ± 1°C
- Temperature stability typically  $\pm 0.1$ °C
- Specifications valid for ambient temperatures from 15°C to 35°C

### 2.6.2.2.5. External Control Possibilities

• USB as standard.

### 2.6.2.2.6. Dimensions

Width: 170 mm Height: 280 mm Depth: 430 mm Weight: 6 kg

### 2.6.2.2.7. Electrical data

- Mains supply: 115/120 V or 220/240 V AC 10% (factory setting 220/240 V)
- Frequency: 50/60 Hz
- Power: 120 VA

### 2.6.2.2.8. Description of Symbols



CAUTION (Refer to accompanying documents).

Symbol identifying the use of substances dangerous for health and environment. For these substances reference must be made to the laws and standards in force in the laboratory where this instrument is used. If using flammable or explosive substances take the necessary precautions.

**Note:** Connection to the protective ground is ensured through the power supply cable which has to be in accordance with CEE 7 (IEC 23-5).

### 2.6.2.2.9. Environmental conditions

Performance and safety of the Controller are guaranteed at all times if it is operated under the following environmental conditions:

- Installation category II with smaller transient over voltages (standard IEC 664)
- Pollution degree 2 (standard IEC EN 61010-1)
- The instrument is designed for internal use and altitudes below 2000 m
- Mains voltage fluctuations must not exceed 10% of the nominal voltage
- In locations where the relative humidity is lower than 80%
- In locations free from dust or vapours of solvents and acids

• Room temperature between 10°C and 35°C

#### 2.6.2.3. User interface

The menu-driven interface of the Controller allows the user to set the temperature and to view the configuration.

Three buttons are used to access the various functions and change the set parameters.

The multi-functional Controller with its display is described below under manual operation.

# 2.6.3. Installation

#### 2.6.3.1. Checking the serial number

The plate containing the serial number is located on the rear panel at the base of the unit. Check that this number is identical to the serial number on the delivery note. If not, please inform Biochrom immediately.

#### 2.6.3.2. Checking mains supply and fuses

The Controller is factory set to operate at 220 VAC.

For operation at a voltage of 120 VAC, the voltage selector at the rear of the

Controller unit must be reset as described below.

Power supply Type of fuse - 250 V

240 VT 1 A, slow-blow

120 VT 2 A, slow-blow

Adjustment for frequency change between 50 and 60 Hz is not necessary.



#### Fig. 3-1: Selecting the voltage

#### 2.6.3.2.1. Changing the Mains Voltage Setting and Fuses

The voltage set is shown in the window of the voltage selector below the mains input connection. To change this setting switch off the unit at the mains switch, remove the mains cable and proceed as follows:

Slide out the fuse holder by inserting a small screwdriver alternatively into the two slots at the left and right and pulling it out of the housing.

Remove the original fuses and replace with those of correct rating (see above) and push the holder back in position.

Move the slider of the voltage selector with a screwdriver to the desired value. The new selected voltage will be now visible.

#### 2.6.3.3. Instrument set-up

#### 2.6.3.3.1. Start up

The Controller is supplied completely assembled. It is only necessary to make the appropriate connections.

Before starting the Controller you should familiarize yourself with the key functions.

- Position the Controller at a suitable position in your system.
- Leave at least 5 cm in front and the back of the Controller for free air circulation.

- Unscrew the tap on the "Vent" connector.
- Mount the "Vent" connector supplied with the instrument.
- Connect the supplied tubes to "Water in" and "Water out" fast connections.
- Connect the tubes to the accessory to be thermostatted.
- Switch on the Controller.
- The instrument initializes and after one minute it is ready to go.
- Check the system for water level ("Water level alarm sensor).

# 2.6.3.3.2. External Connections

The Controller has no special electrical external connections which would allow relay control etc.

Operations are performed either manually or via remote control using the USB connection.

Contact Biochrom for further information about the remote control functions.

# 2.6.4. Operating Instructions

# 2.6.4.1. Manual operation

# Main page and functions of the buttons

Page	ige Display L button Middle button		R button		
Stand	by	Stand by		On	
Main	Normal status	Current temperature Set temperature Heating / cooling power (R)	Menu	Off	Modify the temperature setpoint
Main	Temperature setpoint	Flashing selected parameter	Down	Up	Change
	Alarm status	Alarm description	Menu	Beep on / off	Reset alarm / Change (1)
Menu		Menu selection	Scroll	Select	Exit

(1) Only if the manual alarm reset option has been configured

# Menu page and functions of the buttons

Menu	Description	L button	Middle button	R button
Parameter	User parameters	Scroll	Select	Exit
Configure	Configuration parameters	Scroll	Select	Exit
Diagnostic	Diagnostic values	Scroll	Select	Exit

# Parameters setting

User parameters and functions of the buttons

Number	Description	Default value	L button	Middle button	R button
P-01	Temperature unit of measure	°C	Scroll	Select	Change
P-02	Auto start (after switching on)	NO	Scroll	Select	Change
P-03	Contrast	5	Scroll	Select	Change
P-04	Beep on/off	YES	Scroll	Select	Change
P-05	Manual alarm reset	NO	Scroll	Select	Change

# Configure parameters

A password is required to enter this menu.

CAUTION: all these parameters are correctly set in the factory: no changes should be made by the user.

**Diagnostic parameters** 

Description	Unit of measure
Cell temperature	°C
Radiator temperature	°C
Water resistance	Kohm
Power	%
Cell voltage	Volt

# 2.6.5. Maintenance

The Controller requires very little routine maintenance by the user.

Any operation that requires the units to be opened must only be performed by Biochrom technicians or authorized personnel.

Always unplug from the mains when opening the units.

## 2.6.5.1. Cleaning the instrument

For normal cleaning operations on the instrument only use a cloth dipped in water or neutral detergent, do not use organic or abrasive solvents.

# 2.6.5.2. Water level alarm

Check for signs of water level. It is recommended that this is done prior to starting analysis. If a "Water level" alarm is signalled immediately check for spillage in the accessory compartment and remove it with appropriate measures. Using the refilling bottle supplied to fill the system with liquid via the inlet or outlet connectors until the light turns off.

# 2.6.6. Troubleshooting

The Controller is able to indicate malfunctions and incorrect use. The messages available are directly visible on the display, indicated audibly or transmitted via USB for remote communication.

### 2.6.6.1. Audible alarm

The Controller will send an audible alarm if either a leakage was detected or another failure occurs.

# 2.6.6.1.1. Water level Alarm

The Controller is equipped with a "Water level" sensor. This means even small leakage will be detected.

After the leakage is signalled on the Controller press the RESET ALARM button and remove the leakage immediately by checking all connections and fill up the system with distilled water.

## 2.6.6.1.2. Other system Alarms

The following alarms are available:

- Cell temperature too high
- Radiator temperature too high
- Cell temperature sensor failure
- EEPROM memory failure

If one of these alarms is activated, the Controller will signal the failure by an audible

# signal and display relative messages on the display.

# Action: Switch off the instrument immediately!

If the failure does not disappear after the instrument is switched on again, contact service.

### IMPORTANT!

THE CONTROLLER IS SHIPPED ALREADY FILLED WITH DISTILLED WATER. DURING INSTALLATION, WHEN CONNECTING THE TUBING THE WATER LEVEL IN THE TANK WILL DROP AND WILL NEED TO BE TOPPED UP. ONLY USE THE DISTILLED WATER SUPPLIED IN THE SEPARATE 200 CC CONTAINER INSIDE THE PACKAGING TO TOP UP THE WATER IN THE UNIT.

# 3. ACTIVE ACCESSORIES

Accessories that interact with the instrument's software are named active accessories.

The installation of active accessories that use the internal accessory connector is outlined below.

**Note:** The Ultrospec 7000/8000/9000 will be supplied with a single cell holder as standard; this is removed as described above.

 To insert the accessory, lift the front end, slide into the brackets at the back of the cell chamber (indicated by the two arrows) and tighten the thumb screw by turning clockwise.

**2.** With the accessory inserted and secured using the thumb screw, plug the 15-pin plug into the internal accessory connector.

**3.** With the accessory inserted and plugged into the internal accessory connector, replace front panel and switch on the instrument. The accessory will be calibrated during instrument start-up.

# 3.1. 8-Position Cell Changer 80-7100-08

The 8-position cell changer is designed to speed up analysis by allowing users to automatically measure samples. With a cell changer connected the software will allow the user to perform parallel kinetics measurements.









# 3.2. 5-Position Thermostatted Cell Changer

The 5-position thermostatted cell changer is designed to speed up analysis by allowing users to automatically measure samples that are at a controlled temperature. With a cell changer connected the software will allow the user to perform parallel kinetics measurements.

The 5-position changer is fitted as shown and the water tubing connected to an external water circulator. The inlet from the circulator should be connected to the right hand connector.

## 80-7100-09



# 3.3. Sipper & Flowcell

The Sipper is designed to speed up analysis by allowing users to automatically measure samples by pumping them through a Flowcell using a peristaltic pump head. To install the Sipper carry out the following steps







- Remove the front accessory panel by pulling it upwards
- Plug the Sipper cable into the accessory connector port and slide the Sipper panel into the slot

Biochrom supplies the Sipper with a choice of Tubing kits and Flowcell depending upon customers applications

80-7100-37	Tubing kit for TES Sipper. Includes 6 × Marprene™ pump head tubes and 1 m PTFE waste tubing
80-7100-38	Acid Tubing kit for TES Sipper. Includes 6 × Viton™ pump head tubes and 1 m PTFE waste tubing
80-7100-40	10 mm pathlength Quartz Flowcell 80 μl volume

Spare tubing is available under the following part numbers

80-2080-74	Pump head tubing for Sipper. Includes 6 x Marprene pump head tubes
80-7100-39	Acid pump head tubing for Sipper. Includes 6 x Viton pump head tubes
80-2026-34	PTFE waste tubing
80-2055-13	PTFE tubing with Flowcell connectors

The tubing supplied with the Flowcell from Biochrom is 150 cm in length to allow customers flexibility in how they sip from their sample vessels. Unless you require tubing this long for your application take the tubing with the short Flowcell connector attached and cut the tubing to 65 cm in length.

• Screw the short Flowcell connector into the inlet port on the Flowcell which is marked with an arrow and then feed the tubing through the metal "sample beak" on the Sipper from the inside out.



- Take the tall Flowcell connector and connect it to the outlet port of the Flowcell. Slide the whole Sipper unit up and feed the end of the tube through the hole in the front panel underneath the pump head, once again from the inside out and then slide the Sipper unit back down in it's slot.
- Push the end of this tubing firmly into the end of the supplied pump head tubing. Open the pump head by lifting the front lip and feed in the pump head tubing with the tubing that you have just connected on the left hand side. Close the pump head.



• Take the spare piece of clear tubing with no connectors and push it firmly into the other end of the pump head tube. Put the other end of this into a suitable waste container.

Before using the Sipper to run samples the following checks can be carried out to check it is performing properly – see the relevant software manual for how to do this on your instrument.

- Check for leaks Set the Sipper running in rinse mode, sip water and check that there are no leaks.
- Default air gap the Sipper works by pumping up the selected volume of sample, telling the user to remove the end of the tubing from the sample and the system then pumps again to introduce an air gap between samples. The length of this airgap is set within the software but is dependant upon the length of tubing being used the software default should be OK if you have cut your tubing to 65 cm long, if you are using a different length you may need to change your default air gap.
- Sample volume set up a fixed wavelength method which includes the Sipper and set the sip volume to 1 ml. Fill a 10 ml measuring cylinder, sip several samples and check the volume being sipped is approximately 1 ml, if not recalibrate the Sipper as explained in the relevant software manual.

# 3.3.1. Sipper Troubleshooting

Problems with Sipper systems are generally caused by wrongly set parameters or poor maintenance. Typical faults and their cures include:

# 3.3.1.1. Sample not being sipped

- Check pump head closed
- Check pump head tubing for damage
- Check Flowcell is connected and connectors are tightened

- Check PTFE tubing is pushed firmly into pump head tubing
- Check PTFE tubing is not crushed or damaged
- Check Flowcell for blockages
- Check waste pipe for blockages

# 3.3.1.2. Sample being pumped right through Flowcell

• Check air gap and reset to correct value

## 3.3.1.3. Bubbles in Flowcell

- Check Flowcell connectors are tightened
- Clean Flowcell with detergent solution

## 3.3.1.4. Inconsistent readings

- Check volume being sipped try increasing the volume
- Adjust air gap so the end of the sample is closer to the Flowcell inlet more of the sample flushes the previous sample out of the Flowcell

## 3.3.1.5. Wrong volume being sipped

• Check Sipper calibration

# 4. INSTALLING THE INTERNAL PRINTER

Ensure that the instrument is unplugged before installing the internal printer.

**1.** Before the internal printer can be installed the printer blanking panel must be removed. This is accessed from the back of the unit.

2. Remove the screw securing the printer blanking panel.

**3.** Lift off the printer blanking panel to reveal the connector for the internal printer.

4. Connect the printer to the ribbon cable.









5. Slide the front legs of the printer in place and lower into the correct position.

**6.** Secure to the instrument by screwing into place. Reconnect the instrument to the mains and switch on. Once installed the internal printer will appear in the list





